### **thermo**scientific





# Thermo Scientific Heratherm Heating and Drying Ovens

Efficient, safe and easy



Thermo Fisher S C | E N T | F | C

# Thermo Scientific Heratherm Heating and Drying Ovens

## your samples

We are obsessed with sample integrity. Each of our laboratory heating and drying ovens is designed with sample protection as top priority.



# Heratherm<sup>®</sup> ovens are available in three different models. Choose gravity or mechanical convection technology.

- General Protocol units are optimally suited to everyday heating and drying applications up to 250 °C.
- Advanced Protocol models provide the necessary technologies for users requiring ultimate accuracy and increased heating capability up to 330 °C.
- Advanced Protocol Security ovens offer additional safety features aimed at ultimate protection of your most precious samples and products up to 330 °C.

Sizes include three table top models of 60 L, 100 L and 180 L, and two free standing models of 400 L and 750 L.





# Discover why laboratories across the globe rely on Thermo Scientific products every day.

From simple drying of glassware to the most complex and carefully controlled heating applications, your oven is an invaluable part of your daily work. The Heratherm oven range includes an array of solutions to match your needs – and are backed by a two-year warranty on parts and labor, meaning confidence for your investment.\*

\*Subject to Thermo Fisher Scientific Standard Limited Warranty. See www.thermofisher.com or product literature for details.

# efficient

- Improving energy efficiency, increasing cost effectiveness, reducing environmental impact compared to previous models\*\*
- Unique door design reduces heat emission
- Timer for effective usage and additional energy savings
- Space saving footprint without compromising volume
- Flexible shelving system allows for optimal space utilization

# safe

- Smart safety system surveys functionality of oven consistently
- Temperature uniformity as good as +/-1.3°C at 150°C
- Automatic audible and visual over-temperature alarm
- Access port for data logging with independent sensors
- Low external surface temperature for easy handling

## easy

- Easy-to-read, large vacuum fluorescent display
- Easy setting with intuitive interface
- Easy cleaning with rounded corners

# Innovative technology for outstanding versatility



HERATHERM FEATURES	GENERAL P OVE		ADVANCED F OVE		ADVANCED PROTOCOL SECURITY OVENS
	table top	large capacity	table top	large capacity	table top
Temperature range	50 °C¹ -	250 °C	50 °C¹ - 330 °C {	50¹ °C - 250 °C	50 °C¹ - 330 °C
Interior material	stainless ste	el 1.4016	stainless steel	1.4301	stainless steel 1.4301
Rounded interior corners	V	,	V		<b>✓</b>
Microprocessor control	V	,	V		✓
Automatic overtemperature alarm	V	,	<i>V</i>		✓
Stackable with optional stacking kit	<b>✓</b>	•	<b>✓</b>	•	✓
Easy calibration routine	·	,	<i>V</i>		✓
Gravity convection	choice	<b>✓</b>	choice	•	choice
Mechanical convection	choice	•	choice	<b>✓</b>	choice
Timer functions	on/	off	choice of weekly/ ho	ourly/ real time <sup>2</sup>	choice of weekly/ hourly/ real time <sup>2</sup>
Fresh air damper	manually o	controlled	electronically	controlled	electronically controlled
Access port for independent data monitoring	use exha	ust vent	<i>V</i>		✓
RS232 interface	V	,	<i>V</i>		✓
Adjustable fan speed	•		5 speeds <sup>3</sup>	2 speeds <sup>3</sup>	5 speeds <sup>3</sup>
Programmable (ramps / dwells)	•		<i>V</i>		<b>✓</b>
Boost function for accelerated heat-up	•		<b>✓</b>	•	✓
Optional stainless steel exterior	•		<b>V</b>	<b>✓</b>	✓
Dry alarm contact for connection of alarm dev	vice •		<i>V</i>		<b>✓</b>
Automatic undertemperature alarm	•		•		<b>✓</b>
Door alarm	•		•	<b>V</b>	✓
Lockable door	•		•		<b>✓</b>
Connection for optional sample temperature s	sensor •		•		<b>✓</b>

<sup>&</sup>lt;sup>1</sup> Temperatures as low as ambient +10 °C can be selected – requires open damper and no additional heat in unit. <sup>2</sup> for details on timer see page 8. <sup>3</sup> Fan only on mechanical convection models; fan cannot be switched off, will run at least at ~20%; fan speed depends on temperature - with high temperatures, the fan will run faster to prevent overheating of motor. The fan continues running when opening the door.

#### Convection technologies - your choice

01

#### **Mechanical convection**

- Fast sample drying and heating
- Superior temperature uniformity
- Fast temperature recovery after door opening

02

#### **Gravity convection**

- Gentle drying and heating with low air turbulences
- Perfect for working with materials that are sensitive to airflow such as powders and foils



# Select the right model for your needs

#### Thermo Scientific Heratherm Ovens Selector Guide

APPLICATION	MATERIAL/ SAMPLE*	SAMPLE REQUIREMENTS	RECOMMENDED HERATHERM OVEN
Drying	Glassware	Fast drying	Gravity convection for gentle drying, mechanical convection for
	Powders	Gentle / fast drying	fast drying. Choose Advanced Protocol for highest temperature
	Paper, textile	Gentle / fast drying	<ul> <li>accuracy and selection of fan speed. Choose Advanced Protocol</li> <li>Security for additional sample safety. Check also our Thermo</li> </ul>
	Soil, sand	Gentle drying	Scientific vacuum ovens for speedy drying at low temperatures
Electronics Pharmaceut	Electronics	Gentle drying	www.thermoscientific.com/everylab.
	Pharmaceutical preparations	Gentle drying	
Material testing	Cables	Highest temperature accuracy;	Advanced Protocol models with programming capabilities –
		safe processes	mechanical convection models have highest temperature accuracy.
	Plastics	Highest temperature accuracy;	Advanced Protocol Security models for additional sample safety.
		safe processes	
Curing	Paint	Gentle heating	Gravity convection for gentle heating. Choose Advanced Protocol
	Adhesives	Gentle heating	for high temperature accuracy and Advanced Protocol Security for
	Plastics	Gentle heating	additional sample safety.
	Metals	Gentle heating	Gravity convection for gentle heating, or Advanced Protocol
Heated storage	Pills, drugs	Gentle heating; safe processes	mechanical convection at low fan speed. Choose Advanced     Protocol Security for additional sample safety.
Vulcanization	Rubber	Gentle heating	— I Totocoi occurity for additional sample safety.

#### **Footprint Comparison\***

Making the best use of valuable space in your lab, Heratherm ovens are designed with a very small footprint ratio compared to the interior volume. In addition they are stackable with an optional stacking kit.



For heat treatment and drying processes that involve flammable solvents or potentially explosive atmospheres, please talk to your Thermo Scientific representative.

#### **Maximized space efficiency**



The free standing models enhance flexibility by easily moving around on their casters, which can be locked to ensure stable set-up of the units.

### Thermo Scientific Heratherm General Protocol Ovens

Heratherm General Protocol table top ovens are perfect for routine daily work, providing the ideal heating and drying solution for your applications.

Smooth inner casing with easy to clean rounded corners





# technology for efficiency

- Low energy consumption
- Built in timer
- Excellent footprint/volume ratio
- Three sizes (60 L, 100 L, 180 L)
- Inner chambers made from corrosion-resistant stainless steel with rounded corners for easy cleaning

## convenience for ease of use

- · Large, easy to view vacuum fluorescent display
- Simple, microprocessor-based touch button controls
- Doors can be opened over 180° making these units easy to access
- Automatic overtemperature alarm system to protect samples

#### Specifications table/order numbers General Protocol Ovens

Order number		51028139	51028140	51028141
Model		OGS60	OGS100	OGS180
Convection technology		Gravity convection	Gravity convection	Gravity convection
Temperature range	°C	50 °C¹ - 250 °C	50 °C¹ - 250 °C	50 °C¹ - 250 °C
Spatial temperature deviation	at 150 °C	± 4 °C	± 4.5 °C	±5°C
Temperature deviation over tim	e at 150 °C	± 0.4 °C	± 0.4 °C	± 0.4 °C
Footprint	m² / sqft	0.3 / 3.2	0.36 / 3.9	0.47 / 5.1
Chamber volume	L / cuft	65 / 2.3	105 / 3.75	176 / 6.2
Dimensions chamber	r, mm / in (W x H x D)	328 x 480 x 415 / 12.9 x 18.9 x 16.3	438 x 580 x 414 / 17.2 x 22.8 x 16.3	438 x 680 x 589 / 17.2 x 26.8 x 23.2
exterio	r², mm / in (W x H x D)	530 x 755 x 565 / 20.9 x 29.7 x 22.2	640 x 855 x 565 / 25.2 x 33.7 x 22.2	640 x 955 x 738 / 25.2 x 37.6 x 29.1
Number of shelves	supplied / positions	2 / 13	2 / 16	2 / 19
Max. shelf load	kg / lb	25 / 55	25 / 55	25 / 55
Rated voltage / frequency	V / Hz	230 / 50/60	230 / 50/60	230 / 50/60
Rated power / max. current	W/A	1800 / 7.9	3100 / 13.5	3100 / 13.5
Energy consumption at 150° C	W	194	261	320
Weight	kg / lb	42 / 93	53 / 117	66 / 146

<sup>&</sup>lt;sup>1</sup> Temperatures as low as ambient +10 °C can be selected – requires open damper and no additional heat in unit. <sup>2</sup> Depth does not include handle/display (65mm / 2.6 in.) and distance spacer at rear (80mm / 3.1 in.); height includes the feet (35mm / 1.4 in.) NOTE: All figures in all tables are typical average values for series devices, based on factory standard following norm DIN12880. Please contact us for certification information or IQ/OQ documents.documents.

#### 180° door opening

Allows convenient access to the interior for hassle free cleaning and easy loading of samples.



Standard RS232 interface on all units

Fresh air damper provides control over air exchanges

Operating up to 250 °C



#### performance for proven results

- Operating from 50 °C up to 250 °C
- Gravity convection units are designed to protect delicate samples while offering temperature uniformity of +/- 3.5 °C with temperature stability of +/- 0.4 °C at 150 °C
- Mechanical convection models provide a temperature uniformity of +/-2.9 °C with temperature stability of +/-0.3 °C at 150 °C

#### data logging

- The exhaust vent can be used as an access port for an external temperature sensor
- All Heratherm ovens come standard with a RS232 data interface
- Optional factory installed access ports available

#### Specifications table/order numbers General Protocol Ovens

(	Order number		51028148	51028149	51028150
1	Model		OMS60	OMS100	OMS180
(	Convection technology		Mechanical convection	Mechanical convection	Mechanical convection
-	Temperature range	°C	50 °C¹ - 250 °C	50 °C¹ - 250 °C	50 °C¹ - 250 °C
(	Spatial temperature deviation	at 150 °C	± 3.3 °C	± 3.5 °C	± 2.8 °C
-	Temperature deviation over time	at 150 °C	± 0.3 °C	± 0.4 °C	± 0.3 °C
F	Footprint	m² / sqft	0.3 / 3.2	0.36 / 3.9	0.47 / 5.1
(	Chamber volume	L / cuft	66 / 2.3	104 / 3.65	179 / 6.3
[	Dimensions chamber	, mm / in (W x H x D)	354 x 508 x 368 / 13.9 x 20 x 14.5	464 x 608 x 368 / 18.3 x 23.9 x 14.5	464 x 708 x 543 / 18.3 x 27.9 x 21.4
	exterior <sup>2</sup>	, mm / in (W x H x D)	530 x 755 x 565 / 20.9 x 29.7 x 22.2	640 x 855 x 565 / 25.2 x 33.7 x 22.2	640 x 955 x 738 / 25.2 x 37.6 x 29.1
1	Number of shelves	supplied / positions	2 / 13	2 / 16	2 / 19
1	Max. shelf load	kg / lb	25 / 55	25 / 55	25 / 55
F	Rated voltage / frequency	V / Hz	230 / 50/60	230 / 50/60	230 / 50/60
F	Rated power / max. current	W/A	1400 / 6.1	3060 / 13.3	3060 / 13.3
[	Energy consumption at 150° C	W	316	410	465
١	Weight	kg / lb	42 / 93	53 / 117	66 / 146

<sup>&</sup>lt;sup>1</sup> Temperatures as low as ambient +10 °C can be selected – requires open damper and no additional heat in unit. <sup>2</sup> Depth does not include handle/display (65mm / 2.6 in.) and distance spacer at rear (80mm / 3.1 in.); height includes the feet (35mm / 1.4 in.) NOTE: All figures in all tables are typical average values for series devices, based on factory standard following norm DIN12880. Please contact us for certification information or IQ/OQ documents.

### Thermo Scientific Heratherm Advanced Protocol Ovens

As well as incorporating all the benefits of our General Protocol table top ovens, the Advanced Protocol range boasts additional features providing even more flexibility and accuracy.



100 L model shown

#### maximum efficiency

- Timer extends the automation options available to the user
- > Automate the oven to turn on or off at pre-set times
- > Weekly function allows to set on / off times for each day of the week individually
- > Real time function allows to set timer based on actual time in hours and minutes
- > Timer is perfect to save energy by running the oven just at needed time


#### ultimate performance


- Capable of operating at temperatures up to 330 °C
- Our highest level of temperature uniformity
- Exceptionally low energy consumption 60 L models need just 170 W (gravity convection) and 275 W (mechanical convection) per hour to maintain 150 °C

#### Specifications table/order numbers Advanced Protocol Ovens

Order number (coated exterior	or)	51028142	51028143	51028144
Model		OGH60	OGH100	OGH180
Order number (stainless stee	l exterior)	51028709	51028711	51028712
Model		OGH60 SS	OGH100 SS	OGH180 SS
Convection technology		Gravity convection	Gravity convection	Gravity convection
Temperature range	°C	50 °C¹ - 330 °C	50 °C¹ - 330 °C	50 °C¹ - 330 °C
Spatial temperature deviation	at 150 °C	± 2.5 °C	± 3.0 °C	± 2.5 °C
Temperature deviation over time	at 150 °C	± 0.3 °C	± 0.3 °C	± 0.3 °C
Footprint	m² / sqft	0.3 / 3.2	0.36 / 3.9	0.47 / 5.1
Chamber volume	L / cuft	61 / 2.15	99 / 3.6	168 / 6.05
Dimensions chamber	r, mm / in (W x H x D)	328 x 480 x 389 / 12.9 x 18.9 x 15.3	438 x 580 x 389 / 17.2 x 22.8 x 15.3	438 x 680 x 564 / 17.2 x 26.8 x 22.2
exterior <sup>2</sup>	<sup>2</sup> , mm / in (W x H x D)	530 x 755 x 565 / 20.9 x 29.7 x 22.2	640 x 855 x 565 / 25.2 x 33.7 x 22.2	640 x 955 x 738 / 25.2 x 37.6 x 29.1
Number of shelves	supplied / positions	2 / 13	2 / 16	2 / 19
Max. shelf load	kg / lb	25 / 55	25 / 55	25 / 55
Rated voltage / frequency	V / Hz	230 / 50/60	230 / 50/60	230 / 50/60
Rated power / max. current	W/A	1810 / 7.9	3100 / 13.5	3100 / 13.5
Energy consumption at 150° C	W	170	210	290
Weight	kg / lb	44 / 97	55 / 121	69 / 152
Temperature deviation over time Footprint Chamber volume Dimensions chamber exteriora Number of shelves Max. shelf load Rated voltage / frequency Rated power / max. current Energy consumption at 150° C	at 150 °C  m² / sqft  L / cuft  r, mm / in (W x H x D)  r, mm / in (W x H x D)  supplied / positions  kg / lb  V / Hz  W / A	± 0.3 °C 0.3 / 3.2 61 / 2.15 328 x 480 x 389 / 12.9 x 18.9 x 15.3 530 x 755 x 565 / 20.9 x 29.7 x 22.2 2 / 13 25 / 55 230 / 50/60 1810 / 7.9 170	± 0.3 °C 0.36 / 3.9 99 / 3.6 438 × 580 × 389 / 17.2 × 22.8 × 15.3 640 × 855 × 565 / 25.2 × 33.7 × 22.2 2 / 16 25 / 55 230 / 50/60 3100 / 13.5 210	± 0.3 °C 0.47 / 5.1 168 / 6.05 438 × 680 × 564 / 17.2 × 26.8 × 22 640 × 955 × 738 / 25.2 × 37.6 × 29 2 / 19 25 / 55 230 / 50/60 3100 / 13.5 290

<sup>&</sup>lt;sup>1</sup> Temperatures as low as ambient +10 °C can be selected – requires open damper and no additional heat in unit. <sup>2</sup> Depth does not include handle/display (65mm / 2.6 in.) and distance spacer at rear (80mm / 3.1 in.); height includes the feet (35mm / 1.4 in.) NOTE: All figures in all tables are typical average values for series devices, based on factory standard following norm DIN12880. Please contact us for certification information or IQ/OQ documents.documents.

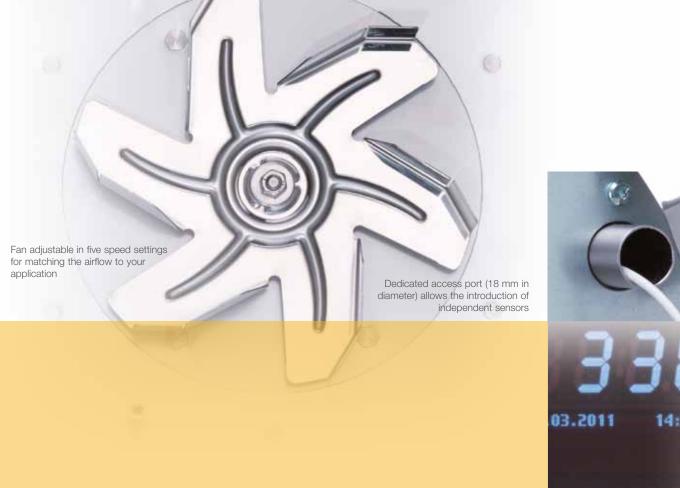

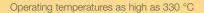



- Adjustable fan speed for application related airflow
- > High speed for fastest drying/heating, quick recovery as well as best temperature stability/uniformity
- > Reduced speed for applications that require minimal airflow, like drying of paper, food or phamaceutical samples
- > Fan speed is adjustable in 5 settings\*
- Programmable controller for temperature ramps and dwells
  - > Store up to 10 programs with 10 discrete steps
  - > Features electronically controlled fan speed and damper position
- Access port allows the introduction of sensors for independent data monitoring
- A simple calibration routine ensures temperature accuracy over time
- Boost function enables rapid heating up no need to run the oven 24/7

\*Actual speed will vary by temperature, and fan cannot be switched off – to prevent overheating

#### Specifications table/order numbers Advanced Protocol Ovens

Order number (coated ext	erior)	51028151	51028152	51028153
Model		OMH60	OMH100	OMH180
Order number (stainless s	teel exterior)	51028713	51028715	51028716
Model		OMH60 SS	OMH100 SS	OMH180 SS
Convection technology		Mechanical convection	Mechanical convection	Mechanical convection
Temperature range	°C	50 °C¹ - 330 °C	50 °C¹ - 330 °C	50 °C1 - 330 °C
Spatial temperature deviation	at 150 °C	± 1.5 °C	± 1.3 °C	± 1.5 °C
Temperature deviation over t	ime at 150 °C	± 0.2 °C	± 0.2 °C	± 0.2 °C
Footprint	m² / sqft	0.3 / 3.2	0.36 / 3.9	0.47 / 5.1
Chamber volume	L / cuft	62 / 2.2	97 / 3.5	170 / 6.1
Dimensions chan	nber, mm / in (W x H x D)	354 x 508 x 343 / 13.9 x 20 x 13.5	464 x 608 x 343 / 18.3 x 23.9 x 13.5	464 x 708 x 518 / 18.3 x 27.9 x 20.4
exte	erior <sup>2</sup> , mm / in (W x H x D)	530 x 755 x 565 / 20.9 x 29.7 x 22.2	640 x 855 x 565 / 25.2 x 33.7 x 22.2	640 x 955 x 738 / 25.2 x 37.6 x 29.1
Number of shelves	supplied / positions	2 / 13	2 / 16	2 / 19
Max. shelf load	kg / lb	25 / 55	25 / 55	25 / 55
Rated voltage / frequency	V / Hz	230 / 50/60	230 / 50/60	230 / 50/60
Rated power / max. current	W/A	1400 / 6.1	3060 / 13.3	3060 / 13.3
Energy consumption at 150°	C W	275	389	448
Weight	kg / lb	44 / 97	55 / 121	69 / 152

<sup>&</sup>lt;sup>1</sup> Temperatures as low as ambient +10 °C can be selected – requires open damper and no additional heat in unit. <sup>2</sup> Depth does not include handle/display (65mm / 2.6 in.) and distance spacer at rear (80mm / 3.1 in.); height includes the feet (35mm / 1.4 in.) NOTE: All figures in all tables are typical average values for series devices, based on factory standard following norm DIN12880. Please contact us for certification information or IQ/OQ documents.

### Thermo Scientific Heratherm Advanced Protocol Security Ovens

Our Advanced Protocol Security table top portfolio combines the benefits of our advanced protocol line with an extra layer of security for applications where process reliability and sample protection are paramount.





All Heratherm Advanced Protocol Security ovens are available in coated or stainless steel exterior. (model 100 L shown)

# enhanced safety

#### Additional security features provide peace-of-mind for precious samples

- Auto-dry function deactivates oven when the samples are dry, saving energy (Note: An optional sample temperature sensor is required to utilize this feature)
- Standard overtemperature alarm and an additional undertemperature alarm ensure that your samples are kept at the correct temperature
- Door lock prevents disruption, tampering or accidental opening
- Door alarm notifies the operator when door is left open accidentally
- Socket for independent sample sensor (option): When connected, exact sample temperature is shown on display for additional peace-of-mind

#### Specifications table/order numbers Advanced Protocol Security Ovens

Order number (coated extended)	erior)	51028145	51028146	51028147
Model		OGH60-S	OGH100-S	OGH180-S
Order number (stainless s	teel exterior)	51028544	51028545	51028546
Model		OGH60-S SS1	OGH100-S SS1	OGH180-S SS <sup>1</sup>
Convection technology		Gravity convection	Gravity convection	Gravity convection
Temperature range	°C	50 °C² - 330 °C	50 °C² - 330 °C	50 °C² - 330 °C
Spatial temperature deviation	at 150 °C	± 2.5 °C	± 2.5 °C	± 2.9 °C
Temperature deviation over t	ime at 150 °C	± 0.3 °C	± 0.3 °C	± 0.3 °C
Footprint	m² / sqft	0.3 / 3.2	0.36 / 3.9	0.47 / 5.1
Chamber volume	L / cuft	61 / 2.15	99 / 3.6	168 / 6.05
Dimensions cham	nber, mm / in (W x H x D)	328 x 480 x 389 / 12.9 x 18.9 x 15.3	438 x 580 x 389 / 17.2 x 22.8 x 15.3	438 x 680 x 564 / 17.2 x 26.8 x 22.2
exte	rior <sup>3</sup> , mm / in (W x H x D)	530 x 755 x 565 / 29.7 x 28.3 x 22.2	640 x 855 x 565 / 25.2 x 33.7 x 22.2	640 x 955 x 738 / 25.2 x 37.6 x 29.1
Number of shelves	supplied / positions	2 / 13	2 / 16	2 / 19
Max. shelf load	kg / lb	25 / 55	25 / 55	25 / 55
Rated voltage / frequency	V / Hz	230 / 50/60	230 / 50/60	230 / 50/60
Rated power / max. current	W/A	1810 / 7.9	3100 / 13.5	3100 / 13.5
Energy consumption at 150°	C W	170	210	290
Weight	kg / lb	44 / 97	55 / 121	69 / 152

¹ Stainless Steel. ² Temperatures as low as ambient +10 °C can be selected – requires open damper and no additional heat in unit. ³ Depth does not include handle/display (65mm / 2.6 in.) and distance spacer at rear (80mm / 3.1 in.); height includes the feet (35mm / 1.4 in.) NOTE: All figures in all tables are typical average values for series devices, based on factory standard following norm DIN12880. Please contact us for certification information or IQ/OQ documents.



An audible door alarm notifies the operator when the door is left open

#### Unique optional sample sensor

- Measure exact sample temperature: shown on display in addition to chamber temperature
- When used with "auto dry" program, the oven will switch off automatically as soon as sample is dry to save energy!





ockable door restricts access

#### Optional stainless steel exterior

- Robust and corrosion-resistant surface
- Easy-to-clean
- Meets demanding needs in pharmaceutical and clinical laboratories



Stainless steel exterior meets demanding needs in pharmaceutical and clinical laboratories

#### Specifications table/order numbers Advanced Protocol Security Ovens

51028154	51028155	51028156
OMH60-S	OMH100-S	OMH180-S
51028547	51028548	51028549
OMH60-S SS1	OMH100-S SS <sup>1</sup>	OMH180-S SS <sup>1</sup>
Mechanical convection	Mechanical convection	Mechanical convection
50 °C² - 330 °C	50 °C² - 330 °C	50 °C² - 330 °C
± 1.5 °C	± 1.3 °C	± 1.5 °C
± 0.2 °C	± 0.2 °C	± 0.2 °C
0.3 / 3.2	0.36 / 3.9	0.47 / 5.1
62 / 2.2	97 / 3.5	170 / 6.1
354 x 508 x 343 / 13.9 x 20 x 13.5	464 x 608 x 343 / 18.3 x 23.9 x 13.5	464 x 708 x 518 / 18.3 x 27.9 x 20.4
530 x 755 x 565 / 20.9 x 29.7 x 22.2	640 x 855 x 565 / 25.2 x 33.7 x 22.2	640 x 955 x 738 / 25.2 x 37.6 x 29.1
2 / 13	2 / 16	2 / 19
25 / 55	25 / 55	25 / 55
230 / 50/60	230 / 50/60	230 / 50/60
1400 / 6.1	3060 / 13.3	3060 / 13.3
275	389	448
44 / 97	55 / 121	69 / 152
	OMH60-S 51028547  OMH60-S SS¹ Mechanical convection 50 °C² - 330 °C ± 1.5 °C ± 0.2 °C 0.3 / 3.2 62 / 2.2 354 x 508 x 343 / 13.9 x 20 x 13.5 530 x 755 x 565 / 20.9 x 29.7 x 22.2 2 / 13 25 / 55 230 / 50/60 1400 / 6.1 275	OMH60-S         OMH100-S           51028547         51028548           OMH60-S SS¹         OMH100-S SS¹           Mechanical convection         Mechanical convection           50 °C² - 330 °C         50 °C² - 330 °C           ± 1.5 °C         ± 1.3 °C           ± 0.2 °C         ± 0.2 °C           0.3 / 3.2         0.36 / 3.9           62 / 2.2         97 / 3.5           354 x 508 x 343 / 13.9 x 20 x 13.5         464 x 608 x 343 / 18.3 x 23.9 x 13.5           530 x 755 x 565 / 20.9 x 29.7 x 22.2         640 x 855 x 565 / 25.2 x 33.7 x 22.2           2 / 13         2 / 16           25 / 55         25 / 55           230 / 50/60         230 / 50/60           1400 / 6.1         3060 / 13.3           275         389

<sup>&</sup>lt;sup>1</sup> Stainless Steel. <sup>2</sup> Temperatures as low as ambient +10 °C can be selected – requires open damper and no additional heat in unit. <sup>3</sup> Depth does not include handle/display (65mm / 2.6 in.) and distance spacer at rear (80mm / 3.1 in.); height includes the feet (35mm / 1.4 in.) NOTE: All figures in all tables are typical average values for series devices, based on factory standard following norm DIN12880. Please contact us for certification information or IQ/OQ documents.

### Thermo Scientific Heratherm Large Capacity General Protocol Ovens

Heratherm large capacity ovens have been designed with your need for larger samples or high sample volume in mind. The General Protocol ovens provide high capacity for day-to-day drying and heating applications.



400 L

750 L

#### efficiency

- Two sizes (400 L and 750 L)
- Gravity convection technology for gentle heating and drying
- Low overall energy consumption: 23% energy savings compared to conventional ovens\*
- Great footprint/volume ratio and optimal use of chamber with flexible shelving system
- Corrosion-resistant stainless steel inner chambers(1.4016)

### exceptional performance

- Operating from 10 °C above ambient temperature up to 250 °C
- Protect delicate samples while still offering a uniformity of +/- 3.0 °C and temperature stability of +/- 0.5 °C at 150 °C

#### ease of use

- "Plug and play" for 230V models with standard plug: no special high voltage connection needed
- Intuitive user interface and easy-to-read display
- Lockable casters for mobility and easy set-up in your lab



Easy-to-view, vacuum fluorescent display with simple-to-use touch button operation

Factory installed casters allow for increased mobility and make setting up your lab easy

Heratherm large capacity ovens offer up to 25%: energy savings compared to conventional ovens\*:





- · Large, easy-to-view, vacuum fluorescent display with simple-to-use touch button operation controlled by an onboard microprocessor
- Doors can be opened to 180° angle for ease-of-access
- Easy-to-remove shelves for cleaning
- Automatic overtemperature alarm system to protect samples no need for timely manual setting
- Timer function to program start or switch off times for additional energy savings
- Lockable casters for easy mobility and stability

#### Specifications table/order numbers Large Capacity General Protocol Ovens

Order number	51029328	51029340	51029342
Model	OGS400	OGS750	OGS750-3P
Convection technology	gravity convection	gravity convection	gravity convection
Temperature range °C	+50 °C¹ - 250 °C	+50 °C¹ - 250 °C	+50 °C¹ - 250 °C
Spatial temperature deviation at 150 °C	±3.0 °C	±3.7 °C	±3.7 °C
Temperature deviation over time at 150 °C	±0.5 °C	±0.5 °C	±0.5 °C
Footprint m² / sqf	0.56 / 6.0	0.91 / 9.8	0.91 / 9.8
Chamber volume L / cuf	419 / 14.8	774 / 27.3	774 / 27.3
Dimensions chamber mm / in (W x H x D	544 x 1307 x 590 / 21.4 x 51.5 x 23.2	1060 x 1335 x 590 / 39.5 x 51.5 x 23.2	1004 x 1307 x 590 / 39.5 x 51.5 x 23.2
exterior <sup>2</sup> mm / in (W x H x D	778 x 1653 x 770 / 30.6 x 66 x 30.3	1261 x 1653 x 770 / 49.6 x 66 x 30.3	1261 x 1653 x 770 / 49.6 x 66 x 30.3
Number of shelves supplied / positions	2/39	2 / 39	2 / 39
Max. shelf load kg / lb	40 / 88	40 / 88	40 / 88
Rated voltage / frequency V / Hz	230 / 50/60	230 / 50/60	400V 3~/N
Rated power / max. current W / A	2420 / 10.5	2990 / 13.0	6350 / 12
Energy consumption at 150° C W	520	800	800
Weight kg / lb	136 / 300	182 / 401	182 / 401

 $<sup>^{1}</sup>$  Temperatures as low as ambient +10  $^{\circ}$ C can be selected – requires open damper and no additional heat in unit

NOTE: All figures in all tables are typical average values for series devices, based on factory standard following norm DIN12880. Please contact us for certification information or IQ/OQ documents.

 $<sup>^2</sup>$  Depth does not include handle/display (65mm / 2.6 in.) and distance spacer at rear (106mm / 4.2 in.); height includes the casters (108mm / 4.3 in.)

### Thermo Scientific Heratherm Large Capacity Advanced Protocol Ovens

The Heratherm Advanced Protocol ovens offer mechanical convection technology for faster drying and better temperature stability and uniformity. Greater flexibility, accuracy and dependability are provided through a wide range of additional features.



# prime performance

- Operating at temperatures as high as 250 °C
- Improved level of temperature uniformity: +/-2.1 to 3.1 °C
- Mechanical convection technology ensures optimal temperature distribution, and fast heat-up and drying processes.
- 2-speed fan for application flexibility:
- > Slow speed for applications that require minimal airflow (e.g. drying of powders),
- > high speed for fastest drying and heating, and best temperature stability and uniformity





#### Additional features

- Access port for introduction of external sensors for independent data monitoring, or other devices for specific test set-ups
- Simple calibration routine to ensure temperature compliance over time.



#### greater efficiency

- Programmable controller for pre-defined temperature ramps
- > Up to 10 temperature steps can be defined per program
- > Fan speed and damper position are controlled electronically, and can be changed for every program step individually for maximal process flexibility
- > Up to 10 programs can be saved for simple and speedy process repetition
- Sophisticated timer extends the automation options available to the user
- > Choose between a simple on/off timer, recurring weekly timer or set oven activity based on the 24 hour clock
- Inner chamber made from stainless steel (highly resistant quality 1.4301)

#### Specifications table/order numbers Large Capacity Advanced Protocol Ovens

Order number (coated exteri	or)	51029331	51029345	51029345
Model		OMH400	OMH750	OMH750-3P
Order number (stainless stee	el exterior)	51029332	51029346	
Model		OMH400-SS	OMH750-SS	n/a
Convection technology		mechanical convection	mechanical convection	mechanical convection
Temperature range	°C	+50 °C¹ - 250 °C	+50 °C¹ - 250 °C²	+50 °C¹ - 300 °C²
Spatial temperature deviation	at 150 °C	±2.1 °C	±3.1 °C	±3.1 °C
Temperature deviation over time	e at 150 °C	±0.3 °C	±0.4 °C	±0.4 °C
Footprint	m² / sqft	0.56 / 6.0	0.91 / 9.8	0.91 / 9.8
Chamber volume	L / cuft	396 / 14	731 / 25.8	731 / 25.8
Dimensions chamber mm	' in (W x H x D)	544 x 1335 x 545 / 21.4 x 52.6 x 21.5	1004 x 1335 x 545 / 39.5 x 52.5 x 21.5	1004 x 1335 x 545 / 39.5 x 52.5 x 21.5
exterior <sup>3</sup> mm	/ in (W x H x D)	778 x 1653 x 770 / 30.6 x 66 x 30.3	1261 x 1653 x 770 / 49.6 x 66 x 30.3	1261 x 1545 x 770 / 49.6 x 60.8 x 30.3
Number of shelves supp	olied / positions	2 / 39	2 / 39	2 / 39
Max. shelf load	kg / lb	40 / 88	40 / 88	40 / 88
Rated voltage / frequency	V / Hz	230 / 50/60	230 / 50/60	400V 3~/N
Rated power / max. current	W/A	2990 / 13.0	3220 / 14.0	6380 / 12.5
Energy consumption at 150° C	W	630	990	990
Weight	kg / lb	135 / 298	185 / 408	185 / 408

<sup>&</sup>lt;sup>1</sup> Temperatures as low as ambient +10 °C can be selected – requires open damper and no additional heat in unit

 $<sup>^2</sup>$  With low voltage and open damper max temperature is 235  $^{\circ}$ C  $^3$  Depth does not include handle/display (65mm / 2.6 in.) and distance spacer at rear (106mm / 4.2 in.); height includes the casters (108mm / 4.3 in.)

### Factory installed options – for table top ovens

#### Specifications table/order numbers

DESCRIPTION	ORDER NUMBER	DETAILS
Options - factory installed: plea	ase order in conjunction	with desired oven
General Protocol kit for	50128887	Kit for underbench installation; for all General Protocol ovens; consists of exhaust tubing
underbench installation		(1.5 m length) and details on wall distances.
Advanced Protocol kit for	51901061	Kit for underbench installation; for all table top Advanced Protocol and Advanced Protocol Security ovens;
underbench installation		consists of exhaust tubing (1.5 m length) and details on wall distances; maximum temperature of oven is set
		at 250 °C for underbench installation. Please order in conjunction with desired oven.
Access port left, small	51900996	Additional access port on center of left side of unit; $\varnothing$ 24 mm for all mechanical convection ovens / $\varnothing$ 19 mm
		for all gravity convection ovens. Please order in conjunction with desired oven.
Access port left, large	51900997	Additional access port on center of left side of unit; $\varnothing$ 58 mm for all mechanical convection ovens / $\varnothing$ 53 mm
		for all gravity convection ovens. Please order in conjunction with desired oven.
Access port right, small	51900998	Additional access port on center of right side of unit; $\varnothing$ 24 mm for all mechanical convection ovens / $\varnothing$ 19 mm
		for all gravity convection ovens. Please order in conjunction with desired oven.
Access port right, large	51900999	Additional access port on center of right side of unit; Ø 58 mm for all mechanical convection ovens / Ø 53 mm
		for all gravity convection ovens. Please order in conjunction with desired oven.
Access port top, small	51901000	Additional access port on center of top of unit; Ø 24 mm for all ovens. Please order in conjunction with
		desired oven.
Access port top, large	51901001	Additional access port on center of top of unit; Ø 58 mm for all ovens. Please order in conjunction with
		desired oven.
Door hinge left side	51900993	Door hinge on left side – available for all table top units. Coated exterior only.
Ovens with view package* - co	ated exterior (stainless	steel exterior not available for this option)
OGH60 with view package	51028897	Advanced Protocol oven with gravity convection; 61 liters volume; 230VAC; 50/60Hz; 1 window; light
OGH100 with view package	51028898	Advanced Protocol oven with gravity convection; 99 liters volume; 230VAC; 50/60Hz; 1 window, light
OGH180 with view package	51028916	Advanced Protocol oven with gravity convection; 168 liters volume; 230VAC; 50/60Hz; 2 windows, light
OMH60 with view package	51028923	Advanced Protocol oven with mechanical convection; 62 liters volume; 230VAC; 50/60Hz; 1 window; light
OMH100 with view package	51028734	Advanced Protocol oven with mechanical convection; 97 liters volume; 230VAC; 50/60Hz; 1 window, light
OMH180 with view package	51028910	Advanced Protocol oven with mechanical convection; 170 I volume; 230VAC; 50/60Hz; 2 windows, light
OGH60-S with view package	51028736	Advanced Protocol Security oven with gravity convection; 61 I volume; 230VAC; 50/60Hz; 1 window; light
OGH100-S with view package	51028924	Advanced Protocol Security oven with gravity convection; 99 I volume; 230VAC; 50/60Hz; 1 window, light
OGH180-S with view package	51028925	Advanced Protocol Security oven with gravity convection; 168 I volume; 230VAC; 50/60Hz; 2 windows, light
OMH60-S with view package	51028926	Advanced Protocol Security oven with mechanical convection; 62 I vol.; 230VAC; 50/60Hz; 1 window; light
OMH100-S with view package	51028927	Advanced Protocol Security oven with mechanical convection; 97 I vol.; 230VAC; 50/60Hz; 1 window, light
OMH180-S with view package	51028733	Advanced Protocol Security oven with mechanical convection; 170 I vol.; 230VAC; 50/60Hz; 2 windows, light

<sup>\*</sup>maximum temperature for ovens with view package is 250 °C

Please ask for our custom solutions. We provide Heratherm table top ovens for cable testing applications, clean room applications, and inner casing with conditional gas-tightness for inert gas applications. Additionally, access ports can be provided for large capacity ovens at preferred positions.



Additional access ports top and right



Underbench installation kit



Viewing windows in 180L unit




Viewing window in 100L unit

### Accessories

#### Specifications table/order numbers

DESCRIPTION ORDER	R NUMBER	DETAILS
Additional shelving - gravity convection		
Additional wire mesh shelf, 60 L models	50127761	For General Protocol; Advanced Protocol; Advanced Protocol Security ovens; incl. 2 shelf supports; 312 x 386 mm
Additional wire mesh shelf, 100 L models	50127762	For General Protocol; Advanced Protocol; Advanced Protocol Security ovens; incl. 2 shelf supports; 422 x 386 mm
Additional wire mesh shelf, 180 L models	50127763	For General Protocol; Advanced Protocol; Advanced Protocol Security ovens; incl. 2 shelf supports; 422 x 561 mm
Additional wire mesh shelf, 400 L models	50135245	For General Protocol ovens; incl. 2 shelf supports; 528 x 535 mm
Additional wire mesh shelf, 750 L models	50135246	For General Protocol ovens; incl. 2 shelf supports; 984 x 535 mm
Reinforced shelf, 180 L models	50128880	Shelf with maximum capacity of 70 kg for Advanced Protocol and Advanced Protocol Security ovens; max total
		capacity is 250 kg
Stainless steel perforated shelf, 400 L model	s 50135241	For General Protocol oven; incl. 2 shelf supports; 528 x 498 mm
Stainless steel perforated shelf, 750 L model	s 50135242	For General Protocol oven; incl. 2 shelf supports; 984 x 498 mm
Additional shelving - mechanical convec	tion	
Additional wire mesh shelf, 60 L models	50127764	For General Protocol; Advanced Protocol; Advanced Protocol Security ovens; incl. 2 shelf supports; 338 x 336 mm
Additional wire mesh shelf, 100 L models	50127765	For General Protocol; Advanced Protocol; Advanced Protocol Security ovens; incl.2 shelf supports; 448 x 336 mm
Additional wire mesh shelf, 180 L models	50127766	For General Protocol; Advanced Protocol; Advanced Protocol Security ovens; incl. 2 shelf supports; 448 x 511 mm
Additional wire mesh shelf, 400 L models	50135245	For Advanced Protocol ovens; incl. 2 shelf supports; 528 x 535 mm
Additional wire mesh shelf, 750 L models	50135246	For Advanced Protocol ovens; incl. 2 shelf supports; 984 x 535 mm
Stainless steel perforated shelf, 60 L models	50127773	For General Protocol, Advanced Protocol, Advanced Protocol Security ovens; incl. 2 shelf supports; 329 x 376 mm
Stainless steel perforated shelf, 100 L mode	ls 50127774	For General Protocol, Advanced Protocol, Advanced Protocol Security ovens; incl. 2 shelf supports; 439 x 376 mm
Stainless steel perforated shelf, 180 L mode	ls 50127777	For General Protocol, Advanced Protocol, Advanced Protocol Security ovens; incl. 2 shelf supports; 439 x 551 mm
Reinforced shelf, 180 L models	50128881	Shelf with maximum capacity of 70 kg for Advanced Protocol and Advanced Protocol Security ovens; max total
		capacity is 250 kg
Stainless steel perforated shelf, 400L models	50135241	For Advanced Protocol oven; incl. 2 shelf supports; 528 x 498 mm
Stainless steel perforated shelf, 750L models	50135242	For Advanced Protocol oven; incl. 2 shelf supports; 984 x 498 mm
Sample Sensor		
Sample sensor, for select models	50127767	Sample sensor for connection to Advanced Protocol Security table top ovens (60 L, 100 L, 180 L); large capacity Advanced Protocol large capacity ovens (400 L, 750 L): measures exact sample temperature; sample temperature is shown on display as plugged in; cable length: 2200 mm
Silicone free viton door sealing		
Silicone free viton door sealing, 60 L models	50130657	Silicone free viton door sealing for all 60 L ovens
Silicone free viton door sealing, 100 L model:	s 50130658	Silicone free viton door sealing for all 100 L ovens
Silicone free viton door sealing, 180 L model:	s 50130659	Silicone free viton door sealing for all 180 L ovens
Silicone free viton door sealing, 400 L model:	s 50135869	Silicone free viton door sealing for all 400 L ovens
Note: there is no silicone-free solution availab	ole for the 750	DL units.
Fresh air filter		
Fresh air particle filter, all ovens models	50127566	Fresh air particle filter for connection to the fresh air vent (intake); for all ovens models
Stacking kits		
Stacking kit, all 60 L models	50126665	Stacking kit to stack two 60 L models
Stacking kit, all 100 L models	50126666	Stacking kit to stack two 100 L models, or 60 L on 100 L
Stacking kit, all 180 L models	50126667	Stacking kit to stack two 100 L models, or 60 L /100 L on 180 L
Support Stands		
Support stand, all 60 L models	50127741	Support stand with casters for 60 L models; height including casters: 187 mm
Support stand, all 100 L models	50127742	Support stand with casters for 100 L models; height including casters: 187 mm
Support stand, all 180 L models	50127743	Support stand with casters for 180 L models; height including casters: 187 mm



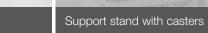





Fresh air particle Filter







LABORATORIUM-APPARATUUR.NL

T. +31 (0)320-266 171 | laboratorium@dijkstra.net pascallaan 9 - 8218 NJ Lelystad | www.dijkstra.net

