

Thermo Scientific CLINIConic™ 30 x 15

Instruction Manual

50121050-d • 07 / 2020

WEEE Conformity

This product is subject to the regulations of the EU Waste Electrical & Electronic Equipment (WEEE) Directive 2012/19/EU. It is marked by the following symbol:

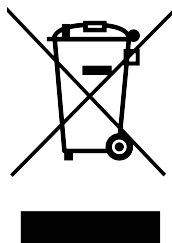


Table of Contents

	Preface	iii
	Scope of Supply	iii
	Precautions	iii
Chapter 1	Rotor Specifications	1-1
	Technical Data	1-2
Chapter 2	Accessories	2-1
Chapter 3	AutoLock™	3-1
	Rotor Installation	3-2
	Removing the Rotor	3-3
Chapter 4	Rotor Loading	4-1
	Before a Run	4-2
	Proper Loading	4-2
	Improper Loading	4-2
	Maximum Loading	4-3
	Cycle Counter	4-4
Chapter 5	Maintenance and Care	5-1
	Cleaning Intervals	5-2
	Cleaning	5-2
	Disinfection	5-3
	Decontamination	5-4
	Autoclaving	5-5
	Service of Thermo Fisher Scientific	5-5
	RCF-Values	A-1
	Chemical Compatibility Chart	B-1

Preface

Before starting to use the rotor, read through the instruction manual carefully and follow the instructions.

The information contained in the instruction manual is the property of Thermo Fisher Scientific; it is forbidden to copy or pass on this information without explicit approval.

Failure to follow the instructions and safety information in this instruction manual will result in the expiration of the seller's warranty.

Scope of Supply

Article Number		Quantity	Check
75003623	CLINIConic™ 30 x 15	1	☐
50121050	Instruction Manual	1	☐
70009824	Anti-corrosion oil	1	☐
76003500	Rubber seal grease	1	☐

If any parts are missing, please contact the nearest Thermo Fisher Scientific representative.

Precautions

In order to ensure safe operation of the CLINIConic™ 30 x 15, the following general safety regulations must be followed:

- Do not remove the magnet at the rotor bottom;
- Do not use rotors which show any signs of corrosion and/or cracks;
- Use only with rotors that have been loaded properly;
- Never run the rotor with missing stainless steel tube holders;
- Never overload the rotor;
- Use only accessories which have been approved by Thermo Fisher Scientific. Exceptions to this rule are commercially available glass or plastic centrifuge tubes, provided they have been approved for the speed or the RCF value of the rotor.

- Please observe the safety instructions.

Please pay particular attention to the following aspects:

- Rotor installation: Check that the rotor is locked properly into place before operating the centrifuge.
- Always balance the samples.

Maximum sample density at maximum speed: $1.2 \frac{\text{g}}{\text{ml}}$



This symbol refers to general hazards.

ATTENTION means that material damage could occur.

WARNING means that injuries or material damage or contamination could occur.



This symbol refers to biological hazards.

Observe the information contained in the instruction manual to keep yourself and your environment safe.

Rotor Specifications

Contents

- “Technical Data” on page 1-2

Technical Data

Table 1-1. 230 V, 50 / 60 Hz ventilated

Centrifuge	Multifuge X1	Megafuge 16	Megafuge 8
Catalog #	75004210	75004230	75007211
Weight empty [kg] of rotor	4.7	4.7	4.7
max. cycle number	50,000	50,000	50,000
Maximum permissible load [g]	30 x 30	30 x 30	30 x 30
maximum speed $n_{\max}$ [rpm]	5,650	5,650	4,400
maximum RCF value at $n_{\max}$	4,997	4,997	3,030
max. / min. radius [cm]	14.0 / 8.5	14.0 / 8.5	14.0 / 8.5
Pitch angle [°]	37	37	37
Accel. / braking time [s]	30 / 35	30 / 35	16 / 27
sample heating at $n_{\max}$ [°C] referred to ambient temperature of 23 °C, running time 60 minutes	9	9	14
aerosol-tight	no	no	no
permissible temperature range autoclavable °C	121	121	121

Table 1-2. 230 V, 50 / 60 Hz refrigerated

Centrifuge	Multifuge X1 R	Megafuge 16R	Megafuge 8R
Catalog #	75004250	75004270	75007214
Weight empty [kg] of rotor	4.7	4.7	4.7
max. cycle number	50,000	50,000	50,000
Maximum permissible load [g]	30 x 30	30 x 30	30 x 30
maximum speed $n_{\max}$ [rpm]	5,650	5,650	4,400
maximum RCF value at $n_{\max}$	4,997	4,997	3,030
max. / min. radius [cm]	14.0 / 8.5	14.0 / 8.5	14.0 / 8.5
Pitch angle [°]	37	37	37
Accel. / braking time [s]	30 / 35	30 / 35	25 / 30
aerosol-tight	no	no	no
permissible temperature range autoclavable °C	121	121	121

Table 1-3. 120 V, 60 Hz ventilated

Centrifuge	Multifuge X1	Megafuge 16	Megafuge 8
Catalog #	75004211	75004231	75007210
Weight empty [kg] of rotor	4.7	4.7	4.7
max. cycle number	50,000	50,000	50,000
Maximum permissible load [g]	30 x 30	30 x 30	30 x 30
maximum speed $n_{\max}$ [rpm]	5,650	5,650	4,400
maximum RCF value at $n_{\max}$	4,997	4,997	3,030
max. / min. radius [cm]	14.0 / 8.5	14.0 / 8.5	14.0 / 8.5
Pitch angle [°]	37	37	37
Accel. / braking time [s]	40 / 40	40 / 40	18 / 26
sample heating at $n_{\max}$ [°C] referred to ambient temperature of 23 °C, running time 60 minutes	9	9	14
aerosol-tight	no	no	no
permissible temperature range autoclavable °C	121	121	121

Table 1-4. 120 V, 60 Hz refrigerated

Centrifuge	Multifuge X1 R	Megafuge 16R	Megafuge 8R
Catalog #	75004251	75004271	75007213
Weight empty [kg] of rotor	4.7	4.7	4.7
max. cycle number	50,000	50,000	50,000
Maximum permissible load [g]	30 x 30	30 x 30	30 x 30
maximum speed $n_{\max}$ [rpm]	5,650	5,650	4,400
maximum RCF value at $n_{\max}$	4,997	4,997	3,030
max. / min. radius [cm]	14.0 / 8.5	14.0 / 8.5	14.0 / 8.5
Pitch angle [°]	37	37	37
Accel. / braking time [s]	40 / 40	40 / 40	25 / 30
aerosol-tight	no	no	no
permissible temperature range autoclavable °C	121	121	121

Table 1-5. 230 V, 50 / 60 Hz ventilated

Centrifuge	Sorvall Legend X1	Sorvall ST 16	Sorvall ST 8
Catalog #	75004220	75004240	75007201
Weight empty [kg] of rotor	4.7	4.7	4.7
max. cycle number	50,000	50,000	50,000
Maximum permissible load [g]	30 x 30	30 x 30	30 x 30
maximum speed $n_{\max}$ [rpm]	5,650	5,650	4,400
maximum RCF value at $n_{\max}$	4,997	4,997	3,030
max. / min. radius [cm]	14.0 / 8.5	14.0 / 8.5	14.0 / 8.5
Pitch angle [°]	37	37	37
Accel. / braking time [s]	30 / 35	30 / 35	16 / 27
sample heating at $n_{\max}$ [°C] referred to ambient temperature of 23 °C, running time 60 minutes	9	9	14
aerosol-tight	no	no	no
permissible temperature range autoclavable °C	121	121	121

Table 1-6. 230 V, 50 / 60 Hz refrigerated

Centrifuge	Sorvall Legend X1R	Sorvall ST 16R	Sorvall ST 8R
Catalog #	75004260	75004380	75007204
Weight empty [kg] of rotor	4.7	4.7	4.7
max. cycle number	50,000	50,000	50,000
Maximum permissible load [g]	30 x 30	30 x 30	30 x 30
maximum speed $n_{\max}$ [rpm]	5,650	5,650	4,400
maximum RCF value at $n_{\max}$	4,997	4,997	3,030
max. / min. radius [cm]	14.0 / 8.5	14.0 / 8.5	14.0 / 8.5
Pitch angle [°]	37	37	37
Accel. / braking time [s]	30 / 35	30 / 35	25 / 30
aerosol-tight	no	no	no
permissible temperature range autoclavable °C	121	121	121

Table 1-7. 120 V, 60 Hz ventilated

Centrifuge	Sorvall Legend X1	Sorvall ST 16	Sorvall ST 8
Catalog #	75004221	75004241	75007200
Weight empty [kg] of rotor	4.7	4.7	4.7
max. cycle number	50,000	50,000	50,000
Maximum permissible load [g]	30 x 30	30 x 30	30 x 30
maximum speed n_{max} [rpm]	5,650	5,650	4,400
maximum RCF value at n_{max}	4,997	4,997	3,030
max. / min. radius [cm]	14.0 / 8.5	14.0 / 8.5	14.0 / 8.5
Pitch angle [°]	37	37	37
Accel. / braking time [s]	40 / 40	40 / 40	18 / 26
sample heating at n_{max} [°C] referred to ambient temperature of 23 °C, running time 60 minutes	9	9	14
aerosol-tight	no	no	no
permissible temperature range autoclavable °C	121	121	121

Table 1-8. 120 V, 60 Hz refrigerated

Centrifuge	Sorvall Legend X1R	Sorvall ST 16R	Sorvall ST 8R
Catalog #	75004261	75004381	75007203
Weight empty [kg] of rotor	4.7	4.7	4.7
max. cycle number	50,000	50,000	50,000
Maximum permissible load [g]	30 x 30	30 x 30	30 x 30
maximum speed n_{max} [rpm]	5,650	5,650	4,400
maximum RCF value at n_{max}	4,997	4,997	3,030
max. / min. radius [cm]	14.0 / 8.5	14.0 / 8.5	14.0 / 8.5
Pitch angle [°]	37	37	37
Accel. / braking time [s]	40 / 40	40 / 40	25 / 30
aerosol-tight	no	no	no
permissible temperature range autoclavable °C	121	121	121

Table 1-9. 100 V, 50 / 60 Hz ventilated

Centrifuge	Sorvall Legend 1X	Sorvall ST 16	Sorvall ST 8
Catalog #	75004223	75004243	75007202
Weight empty [kg] of rotor	4.7	4.7	4.7
max. cycle number	50,000	50,000	50,000
Maximum permissible load [g]	30 x 30	30 x 30	30 x 30
maximum speed $n_{\max}$ [rpm]	5,650	5,650	4,400
maximum RCF value at $n_{\max}$	4,997	4,997	3,030
max. / min. radius [cm]	14.0 / 8.5	14.0 / 8.5	14.0 / 8.5
Pitch angle [°]	37	37	37
Accel. / braking time [s]	45 / 40	45 / 40	18 / 26
sample heating at $n_{\max}$ [°C] referred to ambient temperature of 23 °C, running time 60 minutes	9	9	14
aerosol-tight	no	no	no
permissible temperature range autoclavable °C	121	121	121

Table 1-10. 100 V, 50 / 60 Hz refrigerated

Centrifuge	Sorvall Legend X1R	Sorvall ST 16R	Sorvall ST 8R
Catalog #	75004263	75004383	75007205
Weight empty [kg] of rotor	4.7	4.7	4.7
max. cycle number	50,000	50,000	50,000
Maximum permissible load [g]	30 x 30	30 x 30	30 x 30
maximum speed $n_{\max}$ [rpm]	5,650	5,650	4,400
maximum RCF value at $n_{\max}$	4,997	4,997	3,030
max. / min. radius [cm]	14.0 / 8.5	14.0 / 8.5	14.0 / 8.5
Pitch angle [°]	37	37	37
Accel. / braking time [s]	45 / 40	45 / 40	30 / 30
aerosol-tight	no	no	no
permissible temperature range autoclavable °C	121	121	121

Table 1-11. 230 V, 50 / 60 Hz ventilated

Centrifuge	Thermo Scientific SL 16	Thermo Scientific SL 8
Catalog #	75004000	75007221
Weight empty [kg] of rotor	4.7	4.7
max. cycle number	50,000	50,000
Maximum permissible load [g]	30 x 30	30 x 30
maximum speed $n_{\max}$ [rpm]	5,650	4,400
maximum RCF value at $n_{\max}$	4,997	3,030
max. / min. radius [cm]	14.0 / 8.5	14.0 / 8.5
Pitch angle [°]	37	37
Accel. / braking time [s]	30 / 35	16 / 27
sample heating at $n_{\max}$ [°C] referred to ambient temperature of 23 °C, running time 60 minutes	9	14
aerosol-tight	no	no
permissible temperature range autoclavable °C	121	121

Table 1-12. 230 V, 50 / 60 Hz refrigerated

Centrifuge	Thermo Scientific SL 16R	Thermo Scientific SL 8R
Catalog #	75004030	75007224
Weight empty [kg] of rotor	4.7	4.7
max. cycle number	50,000	50,000
Maximum permissible load [g]	30 x 30	30 x 30
maximum speed $n_{\max}$ [rpm]	5,650	4,400
maximum RCF value at $n_{\max}$	4,997	3,030
max. / min. radius [cm]	14.0 / 8.5	14.0 / 8.5
Pitch angle [°]	37	37
Accel. / braking time [s]	30 / 35	25 / 30
aerosol-tight	no	no
permissible temperature range autoclavable °C	121	121

Table 1-13. 120 V, 60 Hz ventilated

Centrifuge	Thermo Scientific SL 16	Thermo Scientific SL 8
Catalog #	75004001	75007220
Weight empty [kg] of rotor	4.7	4.7
max. cycle number	50,000	50,000
Maximum permissible load [g]	30 x 30	30 x 30
maximum speed $n_{\max}$ [rpm]	5,650	4,400
maximum RCF value at $n_{\max}$	4,997	3,030
max. / min. radius [cm]	14.0 / 8.5	14.0 / 8.5
Pitch angle [°]	37	37
Accel. / braking time [s]	40 / 40	18 / 26
sample heating at $n_{\max}$ [°C] referred to ambient temperature of 23 °C, running time 60 minutes	9	14
aerosol-tight	no	no
permissible temperature range autoclavable °C	121	121

Table 1-14. 120 V, 60 Hz refrigerated

Centrifuge	Thermo Scientific SL 16R	Thermo Scientific SL 8R
Catalog #	75004031	75007223
Weight empty [kg] of rotor	4.7	4.7
max. cycle number	50,000	50,000
Maximum permissible load [g]	30 x 30	30 x 30
maximum speed $n_{\max}$ [rpm]	5,650	4,400
maximum RCF value at $n_{\max}$	4,997	3,030
max. / min. radius [cm]	14.0 / 8.5	14.0 / 8.5
Pitch angle [°]	37	37
Accel. / braking time [s]	40 / 40	25 / 30
aerosol-tight	no	no
permissible temperature range autoclavable °C	121	121

Accessories

Contents

- “Rotor data” on page 2-1
- “Accessories” on page 2-1

CliniConic Fixed Angle Aluminum Rotor

Rotor Specifications

Capacity (ml)	30 x 15 ml
Tube Dimensions (mm)	16,5 x 131
Angle°	37
Maximum Speed (rpm)	5.650
K Factor	3954
Net Weight (kg)	4,7

* Max Speed may vary depending on centrifuge unit.

RCF/Radius	RCF (x g)	Radius (cm)
Maximum	4.997	14
Minimum	3.030	8,5



Cliniconic Fixed Angle Rotor

Cat. No.	Description
75003623	CliniConic Fixed Angle Rotor

Tube Ordering Information

Cat. No.	Tube Vol. (ml)	Vol. Fill (ml)	Description	Qty.	Max. Speed (rpm)	Dims ØxL (mm)	Sealing Assembly Required:			Adapters Required:			Accessories Required:		
							Cat. No.	Qty /Pkg	Description	Cat. No.	Qty /Pkg	No. of Places			
-	15	15	15ml Round / Conical	-	-	16,5 x 131	-	-	-	-	-	-	-	-	-
-	10	-	Blood Collection	-	-	16,5 x 95	-	-	-	75003702	1	1	-	-	-
-	7	-	Blood Collection	-	-	16,5 x 131	-	-	-	11172596	1	1	-	-	-
-	5	-	Blood Collection	-	-	13 x 95	-	-	-	11172595	1	1	-	-	-

AutoLock™

Contents

- “Rotor Installation” on page 3-2
- “Removing the Rotor” on page 3-3

Rotor Installation



CAUTION

Unapproved or incorrectly combined accessories can cause serious damage to the centrifuge.

This rotor is equipped with an AutoLock™-system.

This system is used to automatically lock the rotor to the motor shaft. The rotor does not have to be bolted onto the motor shaft.

Proceed as follows:

1. Open the lid of the centrifuge and if necessary remove any dust, foreign objects or residue from the chamber.
AutoLock™ and O-ring must be clean and undamaged.

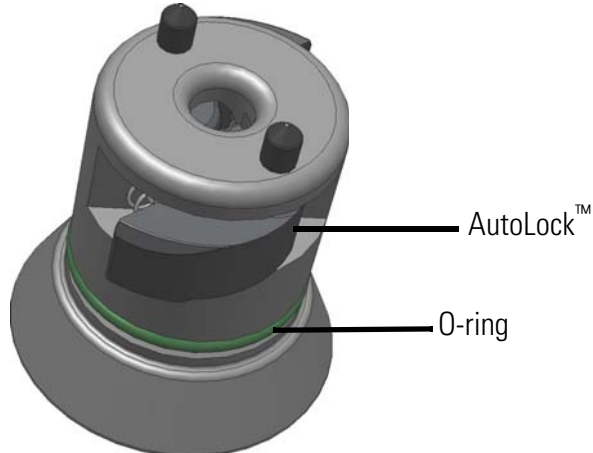


Figure 3-1. AutoLock™

2. Hold the rotor over the shaft and let it slide slowly down the shaft.
The rotor clicks automatically into place.



CAUTION

Do not force the rotor onto the motor shaft.

If the rotor is very light, then it may be necessary to press it onto the shaft with a small amount of pressure.

3. Check if the rotor is properly installed by lifting it slightly on the handle. If the rotor can be pulled up, then it must be reclamped to the shaft.



WARNING

If the rotor cannot be properly locked in place after several attempts, then the AutoLock™ is defective and you are not permitted to operate the rotor.

Check for any damage to the rotor: Damaged rotors must not be used!

Keep the hub area clear of objects.



CAUTION

Check that the rotor is properly locked on the shaft before each use by pulling its handle.



Be sure to check all seals before starting any aerosol-tight applications.
See the information in the rotor instruction manual.

4. Close the centrifuge lid.

Removing the Rotor

To remove the rotor, proceed as follows:

1. Open the centrifuge lid.
2. Grab the rotor handle with both hands and press against the green AutoLock™ key. At the same time, pull the rotor directly upwards with both hands and remove it from the motor shaft. Make sure not to jam the rotor while doing this.

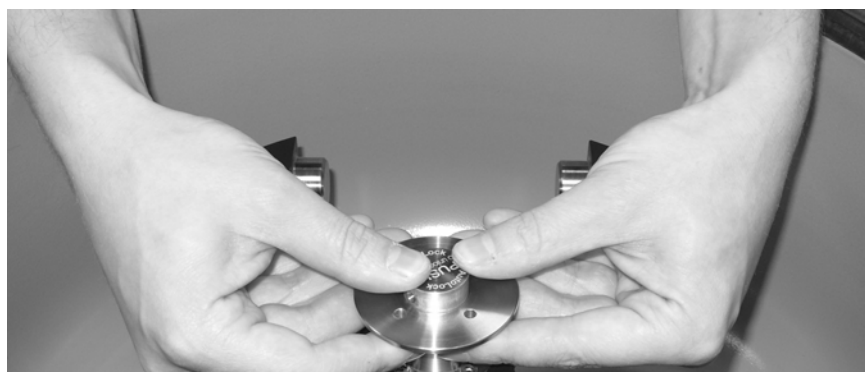


Figure 3-2. Handling the AutoLock™

Rotor Loading

Contents


- “Before a run” on page 4-2
- “Proper Loading” on page 4-2
- “Improper Loading” on page 4-2
- “Maximum loading” on page 4-3
- “Cycle Counter” on page 4-4

Before a Run

1. Please read and observe the safety instructions contained in these operating instructions and in the instructions for use.
2. Check the rotor and all accessory parts for damages such as cracks, scratches or traces of corrosion.
3. Check the rotor chamber, the motor shaft and the AutoLock™.
4. Check the rotor's suitability using the “[Chemical Compatibility Chart](#)” on [page B-1](#).
5. Make sure the tubes or bottles do not touch the lid.


Proper Loading

To ensure safe operation of the centrifuge, the rotor must be evenly loaded at all times.

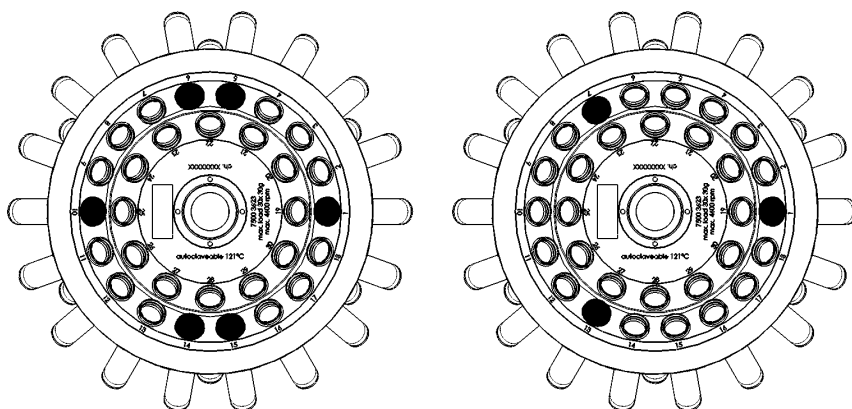

All places must contain stainless steel tube holders. Running the rotor with missing stainless steel tube holders is forbidden. Start loading the rotor on the outer ring. If you only centrifuge one sample place a test tube with water in the opposite position to balance the load.

Improper Loading

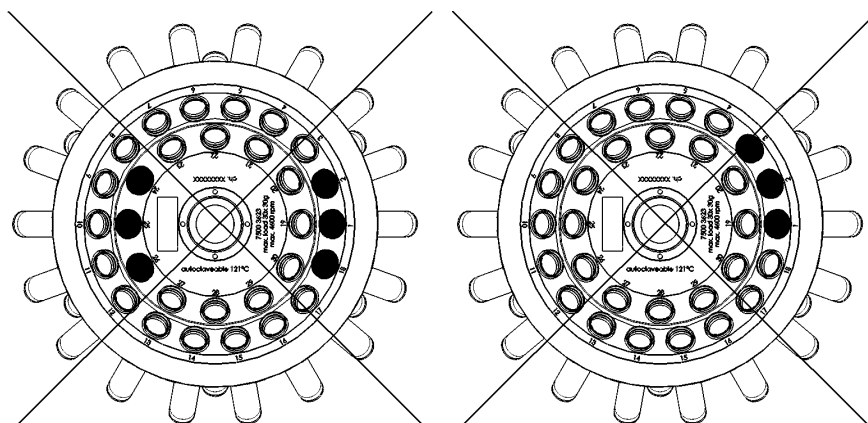

Maximum Loading

The rotor can run at high speeds. The rotor design has sufficient reserve stability even when spinning at top speed.

The safety system of the centrifuge requires that you do not overload the rotor.

There are two options available for centrifuging samples whose weight, including adaptor, exceeds the maximum permissible load:

- Reduce the fill level.
- Reduce the speed.

Use the table or formula below.

Actual Load (g)	RPM _{max}
30	5650
50	4376
70	3699
90	3262
110	2951
130	2714
150	2527
170	2373

Recalculate the maximum permissible speed according to the following formula and select the calculated maximum speed for the centrifuge:

$$n_{\text{adm}} = n_{\text{max}} \sqrt{\frac{\text{Maximum permissible load}}{\text{Effective load}}}$$

n_{adm} = admissible speed

n_{max} = maximum speed

Cycle Counter

The lifetime of rotors is dependent on the amount of mechanical load. Do not exceed the number of cycles recommended for rotors and buckets.

The maximum number of cycles is given in “Rotor Specifications” on [page 1-1](#).



WARNING Replace the rotor when the specified number of cycles is reached. Due to the mechanical load a rotor can break and thus damage the centrifuge.

Service life examples

Usage profile	Maximum service life at 50,000 cycles
frequent use 30 runs / day 220 days / year	7 years

Maintenance and Care

Contents

- “Cleaning Intervals” on page 5-2
- “Cleaning” on page 5-2
- “Disinfection” on page 5-3
- “Decontamination” on page 5-4
- “Autoclaving” on page 5-5
- “Service of Thermo Fisher Scientific” on page 5-5

Cleaning Intervals

For the sake of personal, environmental, and material protection, it is your duty to clean and if necessary, disinfect the centrifuge on a regular basis.

Maintenance	Recommended interval
Clean rotor chamber	daily or when polluted
Clean rotor	daily or when polluted
Accessories	daily or when polluted
Cabinet	Once per month
Ventilation holes	Every six months



CAUTION Refrain from using any other cleaning or decontamination procedure than those recommended here, if you are not entirely sure that the intended procedure is safe for the equipment.
Use only approved cleansers.
If in doubt, contact Thermo Fisher Scientific.

Cleaning

When cleaning centrifuge:

- Use warm water with a neutral solvent.
- Never use caustic cleaning agents such as soap suds, phosphoric acid, bleaching solutions or scrubbing powder.
- Rinse the cavities out thoroughly.
- Use a soft brush without metal bristles to remove stubborn residue.
- Afterwards rinse with distilled water.
- Place the rotors on a plastic grate with their cavities pointing down.
- If drying boxes are used, the temperature must never exceed 50 °C, since higher temperatures could damage the material and shorten the lifetime of the parts.
- Use only disinfectants with a pH of 6-8.
- Dry aluminum parts off with a soft cloth.
- After cleaning, treat the entire surface of aluminum parts with corrosion protection oil (70009824). Also treat the cavities with oil.
- Store the aluminum parts at room temperature or in a cold-storage room with the cavities pointing down.



CAUTION Before using any cleaning or decontamination methods except those recommended by the manufacturer, users should check with the manufacturer that the proposed method will not damage the equipment.

Clean centrifuge and accessories as follows:

1. Open the centrifuge.
2. Turn off the centrifuge.
3. Pull out the power supply plug.
4. Grasp the rotor with both hands and lift it vertically off the centrifuge spindle.
5. Remove the centrifuge tubes and adapters.
6. Use a neutral cleaning agent with a pH value between 6 and 8 for cleaning.
7. Dry all of the rotors and accessories after cleaning with a cloth or in a warm air cabinet at a maximum temperature of 50 °C.
8. Clean the housing of the centrifuge as needed.
 - After cleaning, treat the entire surface of aluminum parts with corrosion protection oil (70009824). Also treat the cavities with oil.
 - Treat the bolt of the swing out rotor with bolt grease (75003786).



CAUTION When cleaning, do not allow liquids, especially organic solvents, to get on the drive shaft, the bearings, the AutoLock™ or the locks. Organic solvents break down the grease in the motor bearing. The drive shaft could freeze up.

After some applications there might be ice in the rotor chamber. Let the ice melt and drain it off. Clean the rotor chamber as described above.

Disinfection

Disinfect the centrifuge immediately whenever infectious material has spilled during centrifugation.



WARNING Infectious material can get into the centrifuge when a tube breaks or as a result of spills. Keep in mind the risk of infection when touching the rotor and take all necessary precautions. In case of contamination, make sure that others are not put at risk. Decontaminate the affected parts immediately. Take other precautions if need be.

Use a sprayer whenever possible so that all surfaces are covered evenly.

The rotor chamber and the rotor should be treated preferably with a neutral disinfectant. A disinfectant spray would be most suitable for this purpose so that the rotor and accessory surfaces are covered evenly.



CAUTION Before using any cleaning or decontamination methods except those recommended by the manufacturer, users should check with the manufacturer that the proposed method will not damage the equipment. Observe the safety precautions and handling instructions for the cleaning agents used.

Contact the Service Department of Thermo Fisher Scientific for questions regarding the use of other disinfectants.

Disinfect the rotor and accessories as follows:

1. Open the centrifuge.
2. Turn off the centrifuge.
3. Pull out the power supply plug.
4. Grasp the rotor with both hands and lift it vertically off the centrifuge spindle.
5. Remove the centrifuge tubes and adapters and dispose of them or disinfect them.
6. Treat the rotor and accessories according to the instructions for the disinfectant (spray or soak in solution). Adhere strictly to the given application times.
7. Be sure the disinfectant can drain off the rotor.
8. Rinse the rotor and rotor lid thoroughly with water and then rub down.
9. Dispose of the disinfectant according to the applicable guidelines.
10. Dry all of the rotors and accessories after cleaning with a cloth or in a warm air cabinet at a maximum temperature of 50 °C.
11. After cleaning, treat the entire surface of aluminum parts with corrosion protection oil (70009824). Also treat the cavities with oil.
12. Treat the bolt of the swing out rotor with bolt grease (75003786).

Decontamination

Decontaminate the centrifuge immediately whenever radioactive material has spilled during centrifugation.



WARNING Radioactive material can get into the centrifuge when a tube breaks or as a result of spills. Keep in mind the risk of infection when touching the rotor and take all necessary precautions.
In case of contamination, make sure that others are not put at risk.
Decontaminate the affected parts immediately.
Take other precautions if need be.



CAUTION Before using any cleaning or decontamination methods except those recommended by the manufacturer, users should check with the manufacturer that the proposed method will not damage the equipment.

For general radioactive decontamination use a solution of equal parts of 70 % ethanol, 10 % SDS and water.

Disinfect the rotor and accessories as follows:

1. Open the centrifuge.
2. Turn off the centrifuge.
3. Pull out the power supply plug.
4. Grasp the rotor with both hands and lift it vertically off the centrifuge spindle.

5. Remove the centrifuge tubes and adapters and dispose of them or disinfect them.
6. Rinse the rotor first with ethanol and then with deionized water.
 - Adhere strictly to the given application times.
7. Be sure the decontamination solution can drain off the rotor.
8. Rinse the rotor and accessories thoroughly with water.
9. Dispose of the decontamination solution according to the applicable guidelines.
10. Dry all of the rotors and accessories after cleaning with a cloth or in a warm air cabinet at a maximum temperature of 50 °C.
11. After cleaning, treat the entire surface of aluminum parts with corrosion protection oil (70009824). Also treat the cavities with oil.
12. Treat the bolt of the swing out rotor with bolt grease (75003786).

Autoclaving

1. Before autoclaving clean rotor and accessories and described above.
2. Place the rotor on a flat surface.
 - Rotors and adapters can be autoclaved at 121 °C.
 - The maximum permissible autoclave cycle is 20 minutes at 121 °C.

Clean the rotor before autoclaving and rinse it with distilled water. Remove all accessories (tubes, adapters) from the rotor. Place the rotor on a flat surface.

Note No chemical additives are permitted in the steam.



CAUTION Never exceed the permitted temperature and duration when autoclaving. If the rotor shows signs of corrosion or wear, it must be replaced.

Service of Thermo Fisher Scientific

Thermo Fisher Scientific recommends having the centrifuge and accessories serviced once a year by an authorized service technician. The service technicians check the following:

- the electrical equipment;
- the suitability of the set-up site;
- the lid lock and the safety system;
- the rotor;
- the fixation of the rotor and the drive shaft.

Thermo Fisher Scientific offers inspection and service contracts for this work.

RCF-Values

Speed (rpm)	R _{min}	R _{max}	RCF R _{min}	RCF R _{max}
300	8.5	14.0	9	14
400	8.5	14.0	15	25
500	8.5	14.0	24	39
600	8.5	14.0	34	56
700	8.5	14.0	47	77
800	8.5	14.0	61	100
900	8.5	14.0	77	127
1000	8.5	14.0	95	157
1100	8.5	14.0	115	189
1200	8.5	14.0	137	225
1300	8.5	14.0	161	265
1400	8.5	14.0	186	307
1500	8.5	14.0	214	352
1600	8.5	14.0	243	401
1700	8.5	14.0	275	452
1800	8.5	14.0	308	507
1900	8.5	14.0	343	565
2000	8.5	14.0	380	626
2100	8.5	14.0	419	690
2200	8.5	14.0	460	758
2300	8.5	14.0	503	828
2400	8.5	14.0	547	902
2500	8.5	14.0	594	978
2600	8.5	14.0	642	1058
2700	8.5	14.0	693	1141
2800	8.5	14.0	745	1227
2900	8.5	14.0	799	1316
3000	8.5	14.0	855	1409

Speed (rpm)	R _{min}	R _{max}	RCF R _{min}	RCF R _{max}
3100	8.5	14.0	913	1504
3200	8.5	14.0	973	1603
3300	8.5	14.0	1035	1705
3400	8.5	14.0	1099	1809
3500	8.5	14.0	1164	1917
3600	8.5	14.0	1232	2028
3700	8.5	14.0	1301	2143
3800	8.5	14.0	1372	2260
3900	8.5	14.0	1445	2381
4000	8.5	14.0	1520	2504
4100	8.5	14.0	1597	2631
4200	8.5	14.0	1676	2761
4300	8.5	14.0	1757	2894
4400	8.5	14.0	1840	3030
4500	8.5	14.0	1924	3170
4600	8.5	14.0	2011	3312
4700	8.5	14.0	2099	3458
4800	8.5	14.0	2189	3606
4900	8.5	14.0	2282	3758
5000	8.5	14.0	2376	3913
5100	8.5	14.0	2472	4071
5200	8.5	14.0	2570	4232
5300	8.5	14.0	2669	4397
5400	8.5	14.0	2771	4564
5500	8.5	14.0	2875	4735
5600	8.5	14.0	2980	4908
5650	8.5	14.0	3034	4997

Chemical Compatibility Chart

CHEMICAL	MATERIAL	ALUMINIUM	ANODIC COATING for ALUMINIUM	BUNA N	CELLULOSE ACETATE BUTYRATE	POLYURETHANE ROTOR PAINT	COMPOSITE Carbon Fiber/Epoxy	DELRIN®	ETHYLENE PROPYLENE	GLASS	NEOPRENE	NORYL®	NYLON	PET®, POLYCLEAR®, CLEARCRIMP®, OCCLEARCRIMP®	POLYALLOMER	POLYCARBONATE	POLYESTER, GLASS THERMOSET	POLYETHERIMIDE	POLYRTHYLENE	POLYPROPYLENE	POLYSULFONE	POLYVINYL CHLORIDE	RULON A®, TEFLON®	SILICONE RUBBER	STAINLESS STEEL	TITANIUM	TYGON®	VITON®
2-mercaptoethanol		S	S	U	-	S	M	S	-	S	U	S	S	U	S	S	-	S	S	S	S	U	S	S	S	S	S	S
Acetaldehyde		S	-	U	U	-	-	-	M	-	U	-	-	-	M	U	U	U	M	M	-	M	S	U	-	S	-	U
Acetone		M	S	U	U	S	U	M	S	S	U	U	S	U	S	U	U	U	S	S	U	U	S	M	M	S	U	U
Acetonitrile		S	S	U	-	S	M	S	-	S	S	U	S	U	M	U	U	-	S	M	U	U	S	S	S	S	U	U
Alconox®		U	U	S	-	S	S	S	-	S	S	S	S	S	S	M	S	S	S	S	S	S	S	S	S	S	S	U
Allyl Alcohol		-	-	-	U	-	-	S	-	-	-	-	S	-	S	S	M	S	S	S	S	-	M	S	-	-	S	-
Aluminum Chloride		U	U	S	S	S	S	U	S	S	S	S	M	S	S	S	S	-	S	S	S	S	S	M	U	U	S	S
Formic Acid (100 %)		-	S	M	U	-	-	U	-	-	-	-	U	-	S	M	U	U	S	S	-	U	S	-	U	S	-	U
Ammonium Acetate		S	S	U	-	S	S	S	-	S	S	S	S	S	S	S	U	-	S	S	S	S	S	S	S	S	S	S
Ammonium Carbonate		M	S	U	S	S	S	S	S	S	S	S	S	S	S	U	U	-	S	S	S	S	S	S	M	S	S	S
Ammonium Hydroxide (10 %)		U	U	S	U	S	S	M	S	S	S	S	S	-	S	U	M	S	S	S	S	S	S	S	S	S	M	S
Ammonium Hydroxide (28 %)		U	U	S	U	S	U	M	S	S	S	S	S	U	S	U	M	S	S	S	S	S	S	S	S	S	M	S
Ammonium Hydroxide (conc.)		U	U	U	U	S	U	M	S	-	S	-	S	U	S	U	U	S	S	S	-	M	S	S	S	S	-	U
Ammonium Phosphate		U	-	S	-	S	S	S	S	S	S	S	S	-	S	S	M	-	S	S	S	S	S	S	M	S	S	S
Ammonium Sulfate		U	M	S	-	S	S	U	S	S	S	S	S	S	S	S	S	-	S	S	S	S	S	S	U	S	S	U
Amyl Alcohol		S	-	M	U	-	-	S	S	-	M	-	S	-	M	S	S	S	S	M	-	-	-	U	-	S	-	M
Aniline		S	S	U	U	S	U	S	M	S	U	U	U	U	U	U	U	-	S	M	U	U	S	S	S	S	U	S
Sodium Hydroxide (<1 %)		U	-	M	S	S	S	-	-	S	M	S	S	-	S	M	M	S	S	S	S	S	S	M	S	S	-	U
Sodium Hydroxide (10 %)		U	-	M	U	-	-	U	-	M	M	S	S	U	S	U	U	S	S	S	S	S	S	M	S	S	-	U
Barium Salts		M	U	S	-	S	S	S	S	S	S	S	S	S	S	S	M	-	S	S	S	S	S	M	S	S	S	S
Benzene		S	S	U	U	S	U	M	U	S	U	U	S	U	U	U	M	U	M	U	U	U	S	U	U	S	U	S
Benzyl Alcohol		S	-	U	U	-	-	M	M	-	M	-	S	U	U	U	U	U	U	U	-	M	S	M	-	S	-	S
Boric Acid		U	S	S	M	S	S	U	S	S	S	S	S	S	S	S	S	U	S	S	S	S	S	S	S	S	S	S

B Chemical Compatibility Chart

CHEMICAL	MATERIAL	POLYMER																														
		ALUMINIUM	ANODIC COATING for ALUMINIUM				BUNA N	CELLULOSE ACETATE BUTYRATE	POLYURETHANE ROTOR PAINT	COMPOSITE Carbon Fiber/Epoxy	DELIRIN®	ETHYLENE PROPYLENE	GLASS	NEOPRENE	NORYL®	NYLON	PET®, POLYCLEAR®, CLEARCRIMP®, CCCLEARCRIMP®	POLYALLUMER	POLYCARBONATE	POLYESTER, GLASS THERMOSET	POLYTHERMIDE	POLYRTHYLENE	POLYPROPYLENE	POLYSULFONE	POLYVINYL CHLORIDE	RULON A®, TEFLON®	SILICONE RUBBER	STAINLESS STEEL	TITANIUM	TYGON®	VITON®	
Cesium Acetate	M	-	S	-	S	S	S	-	S	S	S	S	-	S	S	-	S	S	-	-	S	S	S	S	S	S	S	S	M	S	S	S
Cesium Bromide	M	S	S	-	S	S	S	-	S	S	S	S	S	S	S	S	S	-	-	S	S	S	S	S	S	S	S	M	S	S	S	
Cesium Chloride	M	S	S	U	S	S	S	-	S	S	S	S	S	S	S	S	S	-	-	S	S	S	S	S	S	S	S	M	S	S	S	
Cesium Formate	M	S	S	-	S	S	S	-	S	S	S	S	S	S	S	S	S	-	-	S	S	S	S	S	S	S	S	M	S	S	S	
Cesium Iodide	M	S	S	-	S	S	S	-	S	S	S	S	S	S	S	S	S	-	-	S	S	S	S	S	S	S	S	M	S	S	S	
Cesium Sulfate	M	S	S	-	S	S	S	-	S	S	S	S	S	S	S	S	S	-	-	S	S	S	S	S	S	S	S	M	S	S	S	
Chloroform	U	U	U	U	S	S	M	U	S	U	U	M	U	M	U	U	U	U	U	M	M	U	U	S	U	U	U	M	S			
Chromic Acid (10 %)	U	-	U	U	S	U	U	-	S	S	S	U	S	S	M	U	M	S	S	U	M	S	M	U	S	S	S	S				
Chromic Acid (50 %)	U	-	U	U	-	U	U	-	-	-	S	U	U	S	M	U	M	S	S	U	M	S	-	U	M	-	S					
Cresol Mixture	S	S	U	-	-	-	S	-	S	U	U	U	U	U	U	-	-	U	U	-	U	S	S	S	S	S	U	S				
Cyclohexane	S	S	S	-	S	S	S	U	S	U	S	S	U	U	U	M	S	M	U	M	M	S	U	M	M	U	S					
Deoxycholate	S	S	S	-	S	S	S	-	S	S	S	S	S	S	S	-	-	S	S	S	S	S	S	S	S	S	S	S				
Distilled Water	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S				
Dextran	M	S	S	S	S	S	S	-	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	M	S	S	S		
Diethyl Ether	S	S	U	U	S	S	S	U	S	U	U	S	U	U	U	U	U	U	U	U	U	U	U	S	S	S	S	M	U			
Diethyl Ketone	S	-	U	U	-	-	M	-	S	U	-	S	-	M	U	U	U	M	M	-	U	S	-	-	S	U	U					
Diethylpyrocarbonate	S	S	U	-	S	S	S	-	S	S	U	S	U	S	U	-	-	S	S	S	M	S	S	S	S	S	S	S				
Dimethylsulfoxide	S	S	U	U	S	S	S	-	S	U	S	S	U	S	U	U	-	S	S	U	U	S	S	S	S	U	U					
Dioxane	M	S	U	U	S	S	M	M	S	U	U	S	U	M	U	U	-	M	M	M	U	S	S	S	S	U	U					
Ferric Chloride	U	U	S	-	-	-	M	S	-	M	-	S	-	S	-	-	-	S	S	-	-	-	M	U	S	-	S					
Acetic Acid (Glacial)	S	S	U	U	S	S	U	M	S	U	S	U	U	U	U	U	M	S	U	M	U	S	U	U	S	-	U					
Acetic Acid (5 %)	S	S	M	S	S	S	M	S	S	S	S	S	M	S	S	S	S	S	S	S	S	M	S	S	M	S	S	M				
Acetic Acid (60 %)	S	S	U	U	S	S	U	-	S	M	S	U	U	M	U	S	M	S	M	S	M	S	M	S	M	U	S	M	U			
Ethyl Acetate	M	M	U	U	S	S	M	M	S	S	U	S	U	M	U	U	-	S	S	U	U	S	M	M	S	U	U					
Ethyl Alcohol (50 %)	S	S	S	S	S	S	M	S	S	S	S	S	U	S	U	S	S	S	S	S	S	S	S	S	S	M	S	M	U			
Ethyl Alcohol (95 %)	S	S	S	U	S	S	M	S	S	S	S	S	U	S	U	-	S	S	S	M	S	S	S	U	S	M	U					
Ethylene Dichloride	S	-	U	U	-	-	S	M	-	U	U	S	U	U	U	U	U	U	U	-	U	S	U	-	S	-	S					
Ethylene Glycol	S	S	S	S	S	S	S	S	S	S	S	S	-	S	U	S	S	S	S	S	S	S	S	S	M	S	M	S				
Ethylene Oxide Vapor	S	-	U	-	-	U	-	-	S	U	-	S	-	S	M	-	-	S	S	S	U	S	U	S	S	S	U					
Ficoll-Hypaque®	M	S	S	-	S	S	S	-	S	S	S	S	-	S	S	-	S	S	S	S	S	S	S	M	S	S	S					

CHEMICAL	MATERIAL																													
	ALUMINIUM	ANODIC COATING for ALUMINIUM				BUNA N	CELLULOSE ACETATE BUTYRATE	POLYURETHANE ROTOR PAINT	COMPOSITE Carbon Fiber/Epoxy	DELRI [®]	ETHYLENE PROPYLENE	GLASS	NEOPRENE	NORYL [®]	NYLON	PET ⁺ , POLYCLEAR [®] , CLEARCRIMP [®] , CCCLEARCRIMP [®]	POLYALLUMER	POLYCARBONATE	POLYESTER, GLASS THERMOSET	POLYETHERIMIDE	POLYRTHYLENE	POLYPROPYLENE	POLYSULFONE	POLYVINYL CHLORIDE	RULON A [®] , TEFLON [®]	SILICONE RUBBER	STAINLESS STEEL	TITANIUM	TYGON [®]	VITON [®]
Hydrofluoric Acid (10 %)	U	U	U	M	-	-	U	-	-	U	U	S	-	S	M	U	S	S	S	S	S	M	S	U	U	U	U	-	-	
Hydrofluoric Acid (50 %)	U	U	U	U	-	-	U	-	-	U	U	U	U	S	U	U	U	S	S	M	M	S	U	U	U	U	-	M		
Hydrochloric Acid (conc.)	U	U	U	U	-	U	U	M	-	U	M	U	U	M	U	U	U	-	S	-	U	S	U	U	U	U	-	-		
Formaldehyde (40 %)	M	M	M	S	S	S	S	M	S	S	S	S	M	S	S	S	S	S	S	U	S	S	M	S	S	S	M	S	U	
Glutaraldehyde	S	S	S	S	-	-	S	-	S	S	S	S	S	S	S	S	S	-	-	S	S	S	-	-	S	S	S	-	-	
Glycerol	M	S	S	-	S	S	S	S	S	S	S	S	S	S	S	S	S	S	-	S	S	S	S	S	S	S	S	S	S	
Guanidine Hydrochloride	U	U	S	-	S	S	S	-	S	S	S	S	S	S	S	S	S	-	-	S	S	S	S	S	S	U	S	S	S	
Haemo-Sol [®]	S	S	S	-	-	-	S	-	S	S	S	S	S	S	S	S	S	-	-	S	S	S	S	S	S	S	S	S	S	
Hexane	S	S	S	-	S	S	S	-	S	S	U	S	U	M	U	S	S	U	S	S	U	S	S	M	S	U	S	S	U	
Isobutyl Alcohol	-	-	M	U	-	-	S	S	-	U	-	S	U	S	S	S	M	S	S	S	S	-	S	S	S	-	S	-	S	
Isopropyl Alcohol	M	M	M	U	S	S	S	S	S	U	S	S	U	S	U	S	U	M	S	S	S	S	S	S	S	M	M	M	S	
Iodoacetic Acid	S	S	M	-	S	S	S	-	S	M	S	S	M	S	S	S	-	M	S	S	S	S	S	S	M	S	S	M	M	
Potassium Bromide	U	S	S	-	S	S	S	-	S	S	S	S	S	S	S	S	S	S	S	S	S	S	-	S	S	M	S	S	S	
Potassium Carbonate	M	U	S	S	S	S	S	-	S	S	S	S	S	S	S	U	S	S	S	S	S	S	S	S	S	S	S	S	S	
Potassium Chloride	U	S	S	-	S	S	S	S	S	S	S	S	S	S	S	S	-	S	S	S	S	S	S	S	S	U	S	S	S	
Potassium Hydroxide (5 %)	U	U	S	S	S	S	M	-	S	S	S	S	-	S	U	S	S	S	S	S	S	S	S	S	M	U	M	S	U	
Potassium Hydroxide (conc.)	U	U	M	U	-	-	M	-	M	S	S	-	U	M	U	U	U	S	M	-	M	U	-	U	U	-	U			
Potassium Permanganate	S	S	S	-	S	S	S	-	S	S	S	U	S	S	S	S	M	-	S	M	S	U	S	S	M	S	U	S		
Calcium Chloride	M	U	S	S	S	S	S	S	S	S	S	S	S	S	S	M	S	-	S	S	S	S	S	S	M	S	S	S	S	
Calcium Hypochlorite	M	-	U	-	S	M	M	S	-	M	-	S	-	S	M	S	-	S	S	S	M	S	M	S	M	U	S	-	S	
Kerosene	S	S	S	-	S	S	S	U	S	M	U	S	U	M	M	S	-	M	M	M	S	S	U	S	S	U	S	U	S	
Sodium Chloride (10 %)	S	-	S	S	S	S	S	S	-	-	-	S	S	S	S	S	S	-	S	S	S	S	-	S	S	M	-	S	S	
Sodium Chloride (sat'd)	U	-	S	U	S	S	S	-	-	-	-	S	S	S	S	S	S	-	S	S	-	S	-	S	S	M	-	S	S	
Carbon Tetrachloride	U	U	M	S	S	U	M	U	S	U	U	S	U	M	U	S	S	M	M	S	M	M	M	M	M	U	S	S	S	
Aqua Regia	U	-	U	U	-	-	U	-	-	-	-	-	U	U	U	U	U	U	U	U	U	-	-	-	-	-	S	-	M	
Solution 555 (20 %)	S	S	S	-	-	-	S	-	S	S	S	S	S	S	S	S	-	-	S	S	S	-	S	S	S	S	S	S	S	
Magnesium Chloride	M	S	S	-	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	M	S	S	S	S	
Mercaptoacetic Acid	U	S	U	-	S	M	S	-	S	M	S	U	U	U	U	U	-	S	U	U	S	M	S	U	S	S	S	S	S	
Methyl Alcohol	S	S	S	U	S	S	M	S	S	S	S	S	U	S	U	M	S	S	S	S	S	S	S	S	M	S	M	U	U	
Methylene Chloride	U	U	U	U	M	S	S	U	S	U	U	S	U	U	U	U	U	U	U	M	U	U	U	S	S	M	U	S	U	

B Chemical Compatibility Chart

CHEMICAL	MATERIAL	ALUMINIUM	ANODIC COATING for ALUMINIUM	BUNA N	CELLULOSE ACETATE BUTYRATE	POLYURETHANE ROTOR PAINT	COMPOSITE Carbon Fiber/Epoxy	DELRI [®]	ETHYLENE PROPYLENE	GLASS	NEOPRENE	NORYL [®]	NYLON	PET ⁺ , POLYCLEAR [®] , CLEARCRIMP [®] , CCCLEARCRIMP [®]	POLYALLOMER	POLYCARBONATE	POLYESTER, GLASS THERMOSET	POLYETHERIMIDE	POLYRTHYLENE	POLYPROPYLENE	POLYSULFONE	POLYVINYL CHLORIDE	RULON A [®] , TEFLON [®]	SILICONE RUBBER	STAINLESS STEEL	TITANIUM	TYGON [®]	VITON [®]
Methyl Ethyl Ketone		S	S	U	U	S	S	M	S	S	U	U	S	U	S	U	U	U	S	S	U	U	S	S	S	S	U	U
Metrizamide [®]		M	S	S	-	S	S	S	-	S	S	S	S	-	S	S	-	-	S	S	S	S	S	S	M	S	S	S
Lactic Acid (100 %)		-	-	S	-	-	-	-	-	-	M	S	U	-	S	S	S	M	S	S	-	M	S	M	S	S	-	S
Lactic Acid (20 %)		-	-	S	S	-	-	-	-	-	M	S	M	-	S	S	S	S	S	S	S	M	S	M	S	S	-	S
N-Butyl Alcohol		S	-	S	U	-	-	S	-	-	S	M	-	U	S	M	S	S	S	S	M	M	S	M	-	S	-	S
N-Butyl Phthalate		S	S	U	-	S	S	S	-	S	U	U	S	U	U	U	M	-	U	U	S	U	S	M	M	S	U	S
N, N-Dimethylformamide		S	S	S	U	S	M	S	-	S	S	U	S	U	S	U	U	-	S	S	U	U	S	M	S	S	S	U
Sodium Borate		M	S	S	S	S	S	S	S	S	S	S	U	S	S	S	S	-	S	S	S	S	S	S	M	S	S	S
Sodium Bromide		U	S	S	-	S	S	S	-	S	S	S	S	S	S	S	S	-	S	S	S	S	S	S	M	S	S	S
Sodium Carbonate (2 %)		M	U	S	S	S	S	S	S	S	S	S	S	S	S	U	S	S	S	S	S	S	S	S	S	S	S	S
Sodium Dodecyl Sulfate		S	S	S	-	S	S	S	-	S	S	S	S	S	S	S	-	S	S	S	S	S	S	S	S	S	S	S
Sodium Hypochlorite (5 %)		U	U	M	S	S	M	U	S	S	M	S	S	S	M	S	S	S	S	M	S	S	S	M	U	S	M	S
Sodium Iodide		M	S	S	-	S	S	S	-	S	S	S	S	S	S	S	S	-	-	S	S	S	S	S	M	S	S	S
Sodium Nitrate		S	S	S	-	S	S	S	S	S	S	S	S	S	S	S	S	-	S	S	S	S	S	U	S	S	S	S
Sodium Sulfate		U	S	S	-	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	M	S	S	S
Sodium Sulfide		S	-	S	S	-	-	-	S	-	-	-	S	S	S	U	U	-	-	S	-	-	-	S	S	M	-	S
Sodium Sulfite		S	S	S	-	S	S	S	S	M	S	S	S	S	S	S	M	-	S	S	S	S	S	S	S	S	S	S
Nickel Salts		U	S	S	S	S	S	-	S	S	S	-	-	S	S	S	S	-	S	S	S	S	S	S	M	S	S	S
Oils (Petroleum)		S	S	S	-	-	-	S	U	S	S	S	S	U	U	M	S	M	U	U	S	S	S	U	S	S	S	S
Oils (Other)		S	-	S	-	-	-	S	M	S	S	S	S	U	S	S	S	S	U	S	S	S	S	-	S	S	M	S
Oleic Acid		S	-	U	S	S	S	U	U	S	U	S	S	M	S	S	S	S	S	S	S	S	S	M	U	S	M	M
Oxalic Acid		U	U	M	S	S	S	U	S	S	S	S	S	U	S	U	S	S	S	S	S	S	S	S	U	M	S	S
Perchloric Acid (10 %)		U	-	U	-	S	U	U	-	S	M	M	-	-	M	U	M	S	M	M	-	M	S	U	-	S	-	S
Perchloric Acid (70 %)		U	U	U	-	-	U	U	-	S	U	M	U	U	M	U	U	U	M	M	U	M	S	U	U	S	U	S
Phenol (5 %)		U	S	U	-	S	M	M	-	S	U	M	U	U	S	U	M	S	M	S	U	U	S	U	M	M	M	S
Phenol (50 %)		U	S	U	-	S	U	M	-	S	U	M	U	U	U	U	U	S	U	M	U	U	S	U	U	U	M	S
Phosphoric Acid (10 %)		U	U	M	S	S	S	U	S	S	S	S	U	-	S	S	S	S	S	S	S	S	S	U	M	U	S	S
Phosphoric Acid (conc.)		U	U	M	M	-	-	U	S	-	M	S	U	U	M	M	S	S	S	M	S	M	S	U	M	U	-	S
Physiologic Media (Serum, Urine)		M	S	S	S	-	-	S	-	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S
Picric Acid		S	S	U	-	S	M	S	S	S	M	S	U	S	S	S	U	S	S	S	S	U	S	U	M	S	M	S

CHEMICAL	MATERIAL	ALUMINIUM	ANODIC COATING for ALUMINIUM	BUNA N	CELLULOSE ACETATE BUTYRATE	POLYURETHANE ROTOR PAINT	COMPOSITE Carbon Fiber/Epoxy	DELIN®	ETHYLENE PROPYLENE	GLASS	NEOPRENE	NORYL®	NYLON	PET*, POLYCLEAR®, CLEARCRIMP®, CCCLEARCRIMP®	POLYALLUMER	POLYCARBONATE	POLYESTER, GLASS THERMOSET	POLYTHERMIDE	POLYRTHYLENE	POLYPROPYLENE	POLYSULFONE	POLYVINYL CHLORIDE	RULON A®, TEFLON®	SILICONE RUBBER	STAINLESS STEEL	TITANIUM	TYGON®	VITON®
Pyridine (50 %)		U	S	U	U	S	U	U	-	U	S	S	U	U	M	U	U	-	U	S	M	U	S	S	U	U	U	U
Rubidium Bromide		M	S	S	-	S	S	S	-	S	S	S	S	S	S	S	-	-	S	S	S	S	S	S	M	S	S	S
Rubidium Chloride		M	S	S	-	S	S	S	-	S	S	S	S	S	S	S	-	-	S	S	S	S	S	S	M	S	S	S
Sucrose		M	S	S	-	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S
Sucrose, Alkaline		M	S	S	-	S	S	S	-	S	S	S	S	S	S	U	S	S	S	S	S	S	S	S	M	S	S	S
Sulfosalicylic Acid		U	U	S	S	S	S	S	-	S	S	S	U	S	S	S	-	S	S	S	-	S	S	S	U	S	S	S
Nitric Acid (10 %)		U	S	U	S	S	U	U	-	S	U	S	U	-	S	S	S	S	S	S	S	S	S	M	S	S	S	S
Nitric Acid (50 %)		U	S	U	M	S	U	U	-	S	U	S	U	U	M	M	U	M	M	M	S	S	S	U	S	S	M	S
Nitric Acid (95 %)		U	-	U	U	-	U	U	-	-	U	U	U	U	M	U	U	U	U	M	U	U	S	U	S	S	-	S
Hydrochloric Acid (10 %)		U	U	M	S	S	S	U	-	S	S	S	U	U	S	U	S	S	S	S	S	S	S	S	U	M	S	S
Hydrochloric Acid (50 %)		U	U	U	U	S	U	U	-	S	M	S	U	U	M	U	U	S	S	S	S	M	S	M	U	U	M	M
Sulfuric Acid (10 %)		M	U	U	S	S	U	U	-	S	S	M	U	S	S	S	S	S	S	S	S	S	S	U	U	U	S	S
Sulfuric Acid (50 %)		M	U	U	U	S	U	U	-	S	S	M	U	U	S	U	U	M	S	S	S	S	S	U	U	U	M	S
Sulfuric Acid (conc.)		M	U	U	U	-	U	U	M	-	-	M	U	U	S	U	U	U	M	S	U	M	S	U	U	U	-	S
Stearic Acid		S	-	S	-	-	-	S	M	S	S	S	S	-	S	S	S	S	S	S	S	S	S	M	M	S	S	S
Tetrahydrofuran		S	S	U	U	S	U	U	M	S	U	U	S	U	U	U	-	M	U	U	U	U	S	U	S	S	U	U
Toluene		S	S	U	U	S	S	M	U	S	U	U	S	U	U	U	S	U	M	U	U	U	S	U	S	U	U	M
Trichloroacetic Acid		U	U	U	-	S	S	U	M	S	U	S	U	U	S	M	-	M	S	S	U	U	S	U	U	U	M	U
Trichloroethane		S	-	U	-	-	-	M	U	-	U	-	S	U	U	U	U	U	U	U	U	U	S	U	-	S	-	S
Trichloroethylene		-	-	U	U	-	-	-	U	-	U	-	S	U	U	U	U	U	U	U	U	U	S	U	-	U	-	S
Trisodium Phosphate		-	-	-	S	-	-	M	-	-	-	-	-	-	S	-	-	S	S	S	-	-	S	-	-	S	-	S
Tris Buffer (neutral pH)		U	S	S	S	S	S	S	-	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S
Triton X-100®		S	S	S	-	S	S	S	-	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S
Urea		S	-	U	S	S	S	S	-	-	-	-	S	S	S	M	S	S	S	S	-	S	S	S	M	S	-	S
Hydrogen Peroxide (10 %)		U	U	M	S	S	U	U	-	S	S	S	U	S	S	S	M	U	S	S	S	S	S	S	M	S	U	S
Hydrogen Peroxide (3 %)		S	M	S	S	S	-	S	-	S	S	S	S	S	S	S	S	M	S	S	S	S	S	S	S	S	S	S
Xylene		S	S	U	S	S	S	M	U	S	U	U	U	U	U	U	M	U	M	U	U	U	S	U	M	S	U	S
Zinc Chloride		U	U	S	S	S	S	U	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	U	S	S	S
Zinc Sulfate		U	S	S	-	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S	S
Citric Acid (10 %)		M	S	S	M	S	S	M	S	S	S	S	S	S	S	S	S	M	S	S	S	S	S	S	S	S	S	S

*Polyethyleneterephthalate

Key

- S Satisfactory
- M Moderate attack, may be satisfactory for use in centrifuge depending on length of exposure, speed involved, etc.; suggest testing under actual conditions of use.
- U Unsatisfactory, not recommended.
- Performance unknown; suggest testing, using sample to avoid loss of valuable material.

Chemical resistance data is included only as a guide to product use. Because no organized chemical resistance data exists for materials under the stress of centrifugation, when in doubt we recommend pretesting sample lots.

Index

A	
Accessories	2-1
Aerosol-tight rotors	3-3
Autoclaving	5-5
B	
Before a run	4-2
C	
Care	5-1
Chemical Compatibility Chart	B-1
Cleaning	5-2
Cleaning agents	5-3
Cycle counter	4-4
D	
Decontamination	5-4
Disinfection	5-3
I	
Improper loading	4-2
M	
Maintenance	5-1
Maximum loading	4-3
P	
Precautions	iii
Preface	iii
Proper loading	4-2
R	
RCF-values	A-1
Removing the rotor	3-3
Rotor installation	3-2
S	
Scope of supply	iii
Service technician	5-5

**Thermo Electron LED GmbH**

Zweigniederlassung Osterode
Am Kalkberg, 37520 Osterode am Harz
Germany

thermofisher.com/rotors

© 2009-2020 Thermo Fisher Scientific Inc. All rights reserved.

All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries unless otherwise indicated.

Delrin, TEFLON, and Viton are registered trademarks of DuPont. Noryl is a registered trademark of SABIC. POLYCLEAR is a registered trademark of Hongye CO., Ltd. Hypaque is a registered trademark of Amersham Health As. RULON A and Tygon are registered trademarks of Saint-Gobain Performance Plastics. Alconox is a registered trademark of Alconox. Ficoll is a registered trademark of GE Healthcare. Haemo-Sol is a registered trademark of Haemo-Sol. Triton is a registered trademark of Union Carbide Corporation. Valox is a registered trademark of General Electric Co.

Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.

Shown pictures within the manual are examples and may differ considering the set parameters and language. Pictures of the user interface within the manual are showing the English version as example.

Australia +61 39757 4300

Austria +43 1 801 40 0

Belgium +32 9 272 54 82

China +800 810 5118, +400 650 5118

France +33 2 2803 2180

Germany national toll free
0800 1 536 376

Germany international +49 6184 90 6000

India toll free +1800 22 8374

India +91 22 6716 2200

Italy +39 02 95059 552

Japan +81 3 5826 1616

Korea +82 2 2023 0600

Netherlands +31 76 579 55 55

New Zealand +64 9 980 6700

Nordic / Baltic / CIS countries
+358 10 329 2200

Russia +7 812 703 42 15, +7 495 739 76 41

Singapore +82 2 3420 8700

Spain / Portugal +34 93 223 09 18

Switzerland +41 44 454 12 12

UK / Ireland +44 870 609 9203

USA / Canada +1 866 984 3766

Other Asian Countries +852 3107 7600

Countries not listed +49 6184 90 6000

en

