Bedienungsanleitung
Operating Instructions
Mode d'emploi

- » Super-Star-Exsikkatoren
- » Super-Star-Desiccators
- » Dessiccateur Super-Star

SICCO Super-Star-Exsikkatoren, Artikelnummer: V 1875-07, V 1825-07, V 1935-07, V 1932-01, V 1933-01 und V 1996-08

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OPERATING INSTRUCTIONS


1. Supplied with	13
2. Security Advice	13
3. Taking into operation	14
4. Functionality	17
5. Service / Maintenance	19
6. Disturbances / Repairs	19
7. Waste Disposal	19
8. Specifications	20
9. Accessories	22

SICCO Super-Star-Desiccators

Versions:
Super-Star-Desiccator
Super-Star-Desiccator Vitrum

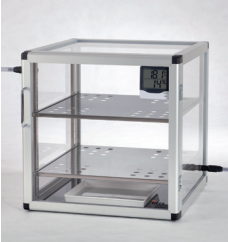SICCO Super Big-Star-Desiccators

Versions:
Super-Big-Star-Desiccator

SICCO Super-Maxi-Desiccators

Versions:
Super-Maxi 1-Desiccator
Super-Maxi 2-Desiccator

SICCO Super Desiccator Wall

Versions:
Super Desiccator Wall

Sicco Super-Star-Desiccators

Please read the information listed below thoroughly. We urgently ask you to observe the instructions about security, use and maintenance.

SICCO Super-Star-Desiccators are made of high-quality and transparent plastic material, except the model Vitrum with panels made of borosilicate glass 3.3. They are made for safe and secure storage of humidity-sensitive products as well as of optical and microelectronic parts in a dust-free atmosphere. This happens by adding SILICA GEL which considerably reduces the air humidity by its absorbing effect or by swapping the SICCO Super-Star-Desiccator with inert gas.

1. Supplied with

	Shelves made of stainless steel	Rails	Hygro-meter	Hose set for gas filling	Desiccant tray	Operating instructions
Super-Star-Desiccator	4	8	1	1	1	1
Super-Star-Desiccator Vitrum	4	8	1	1	1	1
Super-Big-Star-Desiccator	2	4	1	1	1	1
Super-Maxi 1-Desiccator	4	8	1	1	1	1
Super-Maxi 2-Desiccator	4	8	2	2	1	1
Super Desiccator-Wall	6	12	6	0	6	1

2. Security advice

- » Do not take a damaged desiccator into operation.
- » Do not store any solvents or acids in the desiccator except for Super-Star-Desiccator Vitrum.
- » Do not put any explosive or inflammable products into the desiccator.
- » SICCO Super-Star-Desiccators are not suitable for overpressure or vacuum.
- » Provide a safe stand for the desiccator.
- » Place the desiccator in a reachable working height.
- » The ambient temperature range of $-20\text{ }^{\circ}\text{C}$ to $+70\text{ }^{\circ}\text{C}$, respectively of $-20\text{ }^{\circ}\text{C}$ to $+150\text{ }^{\circ}\text{C}$ for Super-Star-Desiccator Vitrum, shall not be exceeded.
- » Do only lead in gas which does not exceed the temperature range of $-20\text{ }^{\circ}\text{C}$ to $+80\text{ }^{\circ}\text{C}$.
- » Do not expose the desiccator to direct sunlight.
- » Remove only one shelf at once.
- » SICCO Super-Star-Desiccators shall be taken into operation only in rooms with sufficient ventilation.
- » For gas filling a connection has to be made to both connectors.
- » Connect the outlet tube at first, then the inlet tube. Disconnection has to be made in reverse order.
- » Make sure that the outlet is open while filling the desiccator with gas to avoid overpressure.
- » Avoid unanticipated reactions which might be caused by gas residues upon opening the door when storing further components in the desiccator.
- » Do not put the filling tubes under tensile or compressive stress.

3. Taking into operation

» Remove the protection cover.

» For models with aluminium rails:
Remove the protection film from the rails.

» Remove the protection film from the shelves.

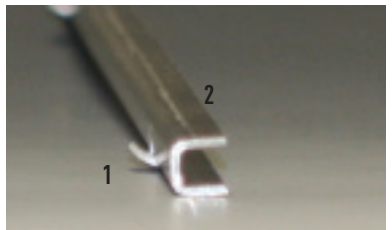
» Opening the door: Hold the desiccator housing with one hand while you open the door by pulling the door grip with the other hand.

» **Please note:** Due to the strong magnetic lock, it is harder to open the door the longer it was closed before. This is normal and can be compensated by pulling stronger than in the everyday use.

» For models with aluminium rails:
Hook the rails (2) into the lateral mounted suspension arrangements (snap-in pin (1) up wards)

» For models with aluminium rails:
Mount the rails (2) horizontally in pairs. Pay attention to mount them in the same height. A slight rotation of the rail helps to fix the snap-in pin safely in the suspension arrangement (3).

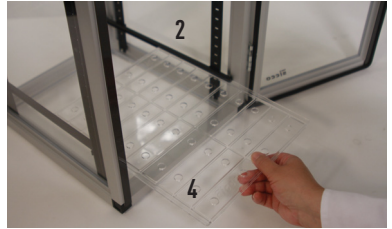
Sicco Super-Star-Desiccators

» For models with rail (2) and suspension arrangement (3) made of plastic:
Mount the rails (2) horizontally in pairs on both sides of the suspension arrangement (3). The engagement of the rail (2) is clearly audible. For a better orientation, the levels are marked by numbers on the suspension arrangement (3).

» The shelves (4) can then be slid in on the rails (2).
» Mount another pair of rails (2) and slide in another shelf (4).
Please note: Mount the rails in pairs and slide in immediately a shelf to avoid that other rails can be fall down accidentally.

» Open the bag with the Silica Gel.
For models Super-Big-Star, Super-Maxi 1 and Super-Maxi 2 the Silica Gel is available as accessorie only.

» Fill the tray (included in delivery) with silica gel. The desiccant an be used immediately.

» Open the battery case on the back of the hygrometer by pushing the cover downwards.
» Remove the battery safety strip and reinsert the battery. Close the battery case.

Assembly of connecting pieces for swapping with gas

» Remove the nut (4) from the connecting piece (6) by turning it counterclockwise. One of the two silicone discs (5) remains on the thread.

Sicco Super-Star-Desiccators

- » One bore is located on the top left, the other one on the bottom right, both in viewing direction to the door. Put the thread of the connecting piece (6) from the outside through the bore. Fix the connecting piece (6) by screwing manually and from the inside the nut (4) on it.

Do not use any tooling for the assembly!

Assemble all other connecting pieces accordingly!

- » Cut the inlet and outlet tubing with a knife to the requested length. 5 m ,respectively 10 m for Super-Maxi 1 and Super-Maxi 2, are included.

- » Press the tubing on the coupling.

Tip: Wet the hose connector with some water. An additional fixation of the tubing is not necessary

- » Pull the slide closure (7) of the coupling (8) in the direction of the tubing and put it on the already mounted connecting piece (6).

Tip: Put some drops of silicone oil onto the sealing surface.

- » For disconnection just hold the coupling (8) and move the slide closure (7) in direction of the coupling.

Attention: The passage on the connecting piece (6) will only be opened after the connection with the coupling has been closed tight. During disconnection, both passages are blocked by an integrated non-return valve.

Please note: gas which is lighter than air has to be led in from the bottom.
Gas which is heavier as air from above. (density of air = 1,2041 kg/m).

Sicco Super-Star-Desiccators

Lockable door (only for Super-Star-Desiccator V 1875-07 and Super-Star-Desiccator Vitrum V 1825-07)

» The lock is open when the cylinder (9) is rebound. Then the key can be removed.

» Putting the cylinder (9) inside and turning the key locks the door. The key can then be removed.

4. Functionality

Rare or inert gases like nitrogen can be introduced as soon as the door is closed tightly. Compared to the ambient air the inert gas does not react with the stored substances.

Alternatively the models Super-Star and Super-Star Vitrum can be dried by using the included silica gel. The silica gel is very hygroscopic and absorbs all air humidity inside the desiccator. The silica gel contains a colour indicator. In unsaturated state, the silica gel is orange. With increasing saturation, the colour is changing from dark brown to black. The silica gel can be regenerated by putting it into an oven for 90 minutes at about 90 to 110 °C. The dumping height influences the time for regeneration: each centimetre in height will take about 90 minutes for regeneration. A microwave oven is not suitable!

Functions of the digital Thermometer/Hygrometer	
Measuring range:	» Temperature – 10 °C to + 60 °C. » Air humidity 20 % to 90 % relative humidity.
Max. error in measurement:	» +/- 1°C at –10 °C to + 50 °C. » +/- 5 % relative humidity at 25 % to 75 % relative humidity and 0 °C to + 50 °C. » Attention: The mounted thermometer/hygrometer cannot display any values or data if the relative inside humidity of the desiccator is less than 25 % or higher than 75 %. To display exact values, we recommend using a thermometer/hygrometer of higher quality.
Batterie:	» For SICCO Star-Desiccator, Star-Desiccator Protect, Star-Desiccator Horizontal, Star-Desiccator Vitrum and Star-Desiccator Antistatic please use a button cell LR44 or AG13. » For SICCO Super-Star-Desiccator please use a battery Micro 1,5 V AAA.

Functionality of the Keys:	Indication of maximum / minimum values: » By pressing the MAX/MIN-key, the maximum value since the last reset will be displayed. » By pressing the MAX/MIN-key again, the minimum value since the last reset will be displayed. » The current values will be displayed after pressing MAX/MIN a third time. Reset of the MAX/MIN-memory: » Press CLEAR and the MAX/MIN-memory will be deleted. The current values will be stored instead. Battery sign in the display » If the battery is low, a battery flag is shown on the left bottom of the display. Exchange the battery to ensure best possible operation of the device. SICCO Super-Star-Desiccator (V1875-07) only: » The key on the back of the thermometer/hygrometer allows to display the temperature either in °C or °F.
-----------------------------------	--

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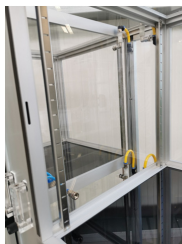
5. Service / Maintenance

- » Use only soft clothes for cleaning. Do not use any detergents which may contain solvents or which scrub.
- » Keep the magnetic door strip clean.

6. Disturbances / Repairs

The following information shall help you to recognise and, if possible, to eliminate possible reasons for disturbances by yourself. If you are not able to handle the disturbance, please give us a call. We will then try to help you this way. If then, a solution to your problem cannot be found, we will ask you to return the cleaned desiccator in the original packing to us. Then we will check whether a cost-effective repair is possible.

Objection	Reason and help
Humidity inside the desiccator does not fall.	» The connecting pieces are not assembled properly. Check again. » The door is not closed tightly. Close the door. » The battery in the hygrometer is empty. Exchange the battery. » The silica gel is saturated. Regenerate the silica gel as described under point 4. » Perhaps you have not added enough silica gel. Add further silica gel. » The built-in hygrometer might be faulty. Double-check the displayed values with an external hygrometer. Exchange the hygrometer.

7. Waste Disposal

The desiccators as well as their packing are made of recyclable materials. Please take further information on the materials used from the following specifications.

8. Specifications

SICCO Super-Star-Desiccators

Cat.-No:	V 1875-07
Material panels:	PMMA (Polymethylmethacrylate)
Temperature resistance:	- 20° to + 70°C
Material frame:	Aluminium with connectors made of polyamide
Material of rails:	Polyamide
Lock:	Magnetic tape
Overall dimensions (W x H x D):	310 x 525 x 375 mm
Additional space required for couplings:	150 mm on each side
Usable interior space (W x H x D):	260 x 480 x 330 mm
Weight:	approx. 7,1 kg
Max. number of shelves:	26 pieces
Max. all-over load per shelf:	10 kg
Total all-over load:	30 kg
Bore dia. of quick couplings:	Ø 6 mm

SICCO Super-Star-Desiccators Vitrum

Cat.-No:	V 1825-07
Material panels:	Borosilicate glass 3.3
Temperature resistance:	- 20° to + 150°C
Material frame:	Aluminium with connectors made of polyamide
Material of rails:	Aluminium
Lock:	Magnetic tape
Overall dimensions (W x H x D):	310 x 525 x 375 mm
Additional space required for couplings:	150 mm on each side
Usable interior space (W x H x D):	260 x 480 x 330 mm
Weight:	approx. 15 kg
Max. number of shelves:	26 pieces
Max. all-over load per shelf:	10 kg
Total all-over load:	30 kg
Bore dia. of quick couplings:	Ø 6 mm

SICCO Super-Big-Star-Desiccators

Cat.-No:	V 1935-07
Material panels:	PMMA (Polymethylmethacrylat)
Temperature resistance:	- 20° to + 70°C
Material frame:	Aluminium with connectors made of polyamide
Material of rails:	Aluminium
Lock:	Magnetic tape
Overall dimensions (W x H x D):	560 x 560 x 580 mm
Additional space required for couplings:	150 mm on each side
Usable interior space (W x H x D):	495 x 500 x 540 mm
Weight:	approx. 18 kg
Max. number of shelves:	17 pieces
Max. all-over load per shelf:	30 kg
Total all-over load:	80 kg
Bore dia. of quick couplings:	Ø 6 mm

SICCO Super-Maxi 1-Desiccators

Cat.-No:	V 1932-07
Material panels:	PMMA (Polymethylmethacrylat)
Temperature resistance:	- 20° to + 70°C
Material frame:	Aluminium with connectors made of polyamide
Material of rails:	Aluminium
Lock:	Magnetic tape
Overall dimensions (W x H x D):	560 x 1150 x 580 mm
Additional space required for couplings:	150 mm on each side
Usable interior space (W x H x D):	495 x 1030 x 540 mm
Weight:	approx. 30 kg
Max. number of shelves:	34 pieces
Max. all-over load per shelf:	30 kg
Total all-over load:	160 kg
Bore dia. of quick couplings:	Ø 6 mm

SICCO Super-Maxi 2-Desiccators

Cat.-No:	V 1933-07
Material panels:	PMMA (Polymethylmethacrylat)
Temperature resistance:	- 20° to + 70°C
Material frame:	Aluminium with connectors made of polyamide
Material of rails:	Aluminium
Lock:	Magnetic tape
Overall dimensions (W x H x D):	560 x 1150 x 580 mm
Additional space required for couplings:	150 mm on each side
Usable interior space (W x H x D):	495 x 500 x 540 mm per chamber
Weight:	approx. 34 kg
Max. number of shelves:	17 pieces per chamber
Max. all-over load per shelf:	30 kg
Total all-over load:	80 kg per chamber
Bore dia. of quick couplings:	Ø 6 mm

SICCO Super-Exsikkatoren Wand

Cat-no.:	V 1996-08
Material panels:	PMMA (Polymethylmethacrylate)
Temperatureresistance:	- 20° to + 70°C
Material frame:	aluminium with connectors made of polyamide
Material of rails:	aluminium
Lock:	magnetictape
Overall dimensions(W x H x D):	1215 x H 1870 x T 580 mm
Usable interior space (W x H x D):	495 x 500 x 540 mm per chamber
Weight:	approx. 86 kg
Max. number of shelves:	17 pc. per chamber
Max. all-over load per shelf.:	30 kg
Total all-over load:	80 kg per chamber
Connection gas supply:	for tubing I.D. 6 mm
Connection exhaustair:	for tubing I.D. Ø 9 mm

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Please note that the Desiccator Wall is not designed for an operation with overpressure or vacuum. Therefore, the exhaust air connection has always to be open during purging with gas.

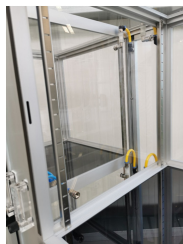
» Before taking into operation close the continuous flow control of each chamber (turn clockwise).

» Then connect the tube for exhaust air to the bigger hose connector and the tube for fresh gas to the smaller connector at the rear side of the desiccator wall. (The tubes are not included in the scope of delivery).

» If one or more compartments are to be filled with gas, open the external gas supply. Then turn the flow meter of the corresponding compartment counter clockwise until the scale shows the required litres per minute. By turning it in clockwise direction, the flow meter is closed again.

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BOHLENDER GmbH