

THE SECRET CODE

OF VACUUM TECHNOLOGY

.....
Good user practice for diaphragm pump systems


MAXIMIZE SERVICE LIFE AND HASSLE-FREE OPERATION



PROTECT THE PUMP MECHANICS BY PREVENTING PARTICLE ASPIRATION
+ Inlet separator



KEEP THE PUMP IN GOOD PERFORMANCE
+ Gas ballast purges condensate and vapors

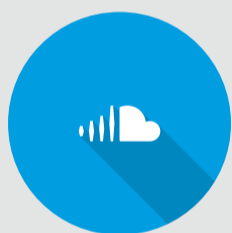

NEVER BLOCK THE PUMPS' OUTLET
+ Free pump exhaust



KEEP THE PUMP CLEAN
+ Post-operation with open gas ballast for "self-cleaning"



MINIMIZE CHEMICAL EMISSION FOR USER HEALTH AND ENVIRONMENTAL PROTECTION
+ Emission condenser for solvent recovery



FOR HASSLE-FREE OPERATION
+ Regular checking and cleaning
+ Maintenance to restore original specs